

# User's Manual

## SteamScrubber<sup>®</sup> & FlaskScrubber<sup>®</sup> Glassware Washers

### Models

|              |              |
|--------------|--------------|
| 44003 Series | 44203 Series |
| 44004 Series | 44204 Series |

To receive important product updates,  
complete your product registration card  
online at **register.labconco.com**



*Protecting your  
laboratory environment*

**LABCONCO**<sup>®</sup>

Labconco Corporation  
8811 Prospect Avenue  
Kansas City, MO 64132-2696  
800-821-5525, 816-333-8811  
FAX 816-363-0130  
E-MAIL [labconco@labconco.com](mailto:labconco@labconco.com)  
HOME PAGE [www.labconco.com](http://www.labconco.com)

**Please read the User's Manual before operating the equipment.**

The information contained in this manual and the accompanying products are copyrighted and all rights reserved by Labconco Corporation. Labconco Corporation reserves the right to make periodic design changes without obligation to notify any person or entity of such change.

## **Warranty**

Labconco provides a warranty on all parts and factory workmanship. The warranty includes areas of defective material and workmanship, provided such defect results from normal and proper use of the equipment.

The warranty for all Labconco products will expire one year from date of installation or two years from date of shipment from Labconco, whichever is sooner, except the following;

- Purifier® Delta® Series Biological Safety Cabinets and PuriCare® Lab Animal Research Stations carry a three-year warranty from date of installation or four years from date of shipment from Labconco, whichever is sooner.
- SteamScrubber® & FlaskScrubber® Glassware Washers carry a two-year warranty from date of installation or three years from date of shipment from Labconco, whichever is sooner.
- Blood Drawing Chairs carry a ten year warranty.
- Carts carry a lifetime warranty.
- Glassware is not warranted from breakage when dropped or mishandled.

This limited warranty covers parts and labor, but not transportation and insurance charges. In the event of a warranty claim, contact Labconco Corporation or the dealer who sold you the product. If the cause is determined to be a manufacturing fault, the dealer or Labconco Corporation will repair or replace all defective parts to restore the unit to operation. Under no circumstances shall Labconco Corporation be liable for indirect, consequential, or special damages of any kind. This statement may be altered by a specific published amendment. No individual has authorization to alter the provisions of this warranty policy or its amendments. Lamps and filters are not covered by this warranty. Damage due to corrosion or accidental breakage is not covered.

## **Returned or Damaged Goods**

Do not return goods without the prior authorization from Labconco. Unauthorized returns will not be accepted. If your shipment was damaged in transit, you must file a claim directly with the freight carrier. Labconco Corporation and its dealers are not responsible for shipping damages.

The United States Interstate Commerce Commission rules require that claims be filed with the delivery carrier within fifteen (15) days of delivery.

## **Limitation of Liability**

The disposal and/or emission of substances used in connection with this equipment may be governed by various federal, state, or local regulations. All users of this equipment are required to become familiar with any regulations that apply in the user's area concerning the dumping of waste materials in or upon water, land, or air and to comply with such regulations. Labconco Corporation is held harmless with respect to user's compliance with such regulations.

## **Contacting Labconco Corporation**

If you have questions that are not addressed in this manual, or if you need technical assistance, contact Labconco's Customer Service Department or Labconco's Product Service Department at 1-800-821-5525 or 1-816-333-8811, between the hours of 7:00 a.m. and 6:00 p.m., Central Standard Time.

---

# TABLE OF CONTENTS

|  |    |
|--|----|
| <b>CHAPTER 1: INTRODUCTION</b>           | 1  |
| Safety Symbols                           | 1  |
| <b>CHAPTER 2: PREREQUISITES</b>          | 3  |
| Hot Water Requirements                   | 3  |
| Purified Water Requirements              | 3  |
| Electrical Requirements                  | 4  |
| Drain Requirements                       | 5  |
| Space Requirements                       | 5  |
| <b>CHAPTER 3: GETTING STARTED</b>        | 6  |
| Unpacking Your Glassware Washer          | 7  |
| Washer Components                        | 7  |
| Removing the Shipping Skid               | 9  |
| Washer Installation                      | 9  |
| Installing the Leveling Feet             | 9  |
| Leveling the Washer                      | 10 |
| Centering the Washer Door                | 11 |
| Securing the Washer to the Countertop    | 11 |
| Connecting the Water Services            | 12 |
| Connecting the Hot Water Supply          | 12 |
| Connecting the Purified Water Supply     | 13 |
| Connecting the Drain Hose                | 15 |
| Connecting the Power Supply              | 17 |
| Controller Set Up                        | 18 |
| Testing Your Washer                      | 19 |
| Running the Diagnostics                  | 20 |
| <b>CHAPTER 4: USING YOUR WASHER</b>      | 22 |
| Arranging the Washer Racks               | 22 |
| The SteamScrubber Racks                  | 22 |
| Removing the Upper Rack                  | 23 |
| The FlaskScrubber Spindle Rack           | 23 |
| Loading Glassware into the FlaskScrubber | 24 |

|  |           |
|--|-----------|
| The Detergent Dispenser                              | 25        |
| Filling the Neutralizing Acid Dispenser              | 26        |
| Selecting an Operating Cycle                         | 27        |
| Programming and Starting a Wash Cycle                | 27        |
| Interrupting a Wash Cycle                            | 29        |
| Alarms   | 30        |
| <b>CHAPTER 5: MAINTAINING YOUR WASHER</b>            | <b>31</b> |
| Maintaining the Stainless Steel Interior             | 31        |
| Maintaining the Air Gap in Drain                     | 32        |
| Maintaining the Water Fill Valve                     | 32        |
| Replacing the Washer Light Bulb                      | 33        |
| Removing the Washer Top, Back, and Side Panels       | 34        |
| Cleaning the Sump Filter Screen                      | 35        |
| <b>CHAPTER 6: MODIFYING YOUR WASHER</b>              | <b>36</b> |
| Adjusting the Height of the Upper Rack               | 36        |
| Spindle Rack Options                                 | 37        |
| Installing Optional Pipet Inserts in a FlaskScrubber | 37        |
| Installing the Optional Upper Spindle Rack           | 38        |
| Installing Optional Inserts into the Spindle Rack    | 40        |
| Installing Small Spindles                            | 41        |
| Installing the Optional Cool Drain Water Kit         | 41        |
| Installing the Optional Liquid Detergent Dispenser   | 42        |
| <b>CHAPTER 7: TROUBLESHOOTING</b>                    | <b>45</b> |
| <b>APPENDIX A: GLASSWARE WASHER ACCESSORIES</b>      | <b>52</b> |
| FlaskScrubber Accessories                            | 52        |
| FlaskScrubber or SteamScrubber Accessories           | 54        |
| Consumables  | 59        |
| <b>APPENDIX B: GLASSWARE WASHER DIMENSIONS</b>       | <b>60</b> |
| <b>APPENDIX C: GLASSWARE WASHER SPECIFICATIONS</b>   | <b>62</b> |
| <b>DECLARATION OF CONFORMITY</b>                     | <b>69</b> |

---

# CHAPTER 1

## INTRODUCTION


Congratulations on your purchase of a Labconco Glassware Washer. Labconco manufactures two types of Glassware Washers, the SteamScrubber® and the FlaskScrubber®. Each of these washers is available in an Undercounter and Freestanding model. An optional kit is available to convert Freestanding models to Mobile models. Each washer model is available in 115V or 230V, and with or without a window in the door.

Your Labconco Glassware Washer is designed and manufactured to thoroughly clean your laboratory glassware and accessories and complies with NSF/ANSI 3 standards for sanitization. You can select pre-programmed wash cycles or program customized wash cycles to meet your requirements.


The Glassware Washer graphic display provides operational information. Models with a window and light in the door allow you to visually monitor the operation of your washer. Wash cycles may be programmed to provide steam and the use of purified water in the rinse cycles, allowing you to choose the best conditions for cleaning your glassware.




### Safety Symbols

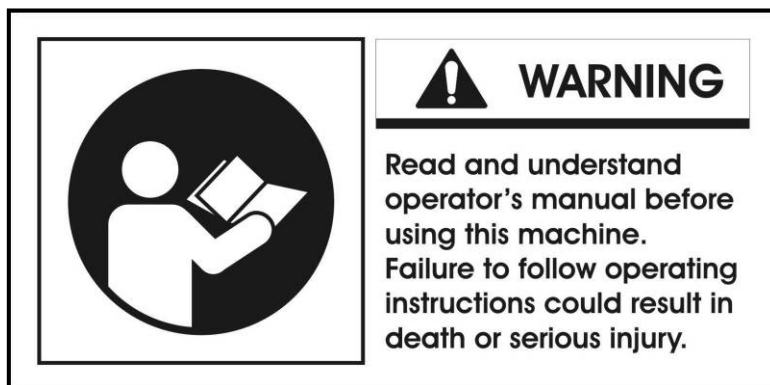Your Glassware Washer was designed with safety in mind, however conditions may exist that could be hazardous.



Throughout this manual potentially hazardous conditions are identified using the following words and symbols.

|   |   |
|---|---|
|  | <p>This is the safety alert symbol. It is used to alert you to potential personal injury hazards. Obey all safety messages that follow this symbol to avoid possible injury or death.</p> |
|---|---|

|  |  |
|--|--|
|  <b>DANGER</b>  | <p><b>DANGER</b> indicates an imminently hazardous situation which, if not avoided, will result in death or serious injury.</p>                              |
|  <b>WARNING</b> | <p><b>WARNING</b> indicates a potentially hazardous situation which, if not avoided, could result in death or serious injury.</p>                            |
|  <b>CAUTION</b> | <p><b>CAUTION</b> indicates a potentially hazardous situation which, if not avoided, may result in minor or moderate injury.</p>                             |
| <b>CAUTION</b>   | <p><b>CAUTION</b> used without the safety alert symbol indicates a potentially hazardous situation which, if not avoided, may result in property damage.</p> |

It is important that you understand the warnings listed throughout this manual before you operate the Glassware Washer.

---

# CHAPTER 2

## PREREQUISITES

Before you install the Glassware Washer, prepare your site for installation. Carefully examine the location where you intend to install the Glassware Washer. You must be certain that the area is level and of solid construction. In addition, a hot water source, a drain, a purified water source (if applicable), and an electrical source must be located near the installation site. A cold water source is required when the optional Cool Drain Water Kit is installed.

Refer to *Appendix C: Glassware Washer Specifications* for complete Glassware Washer electrical and environmental conditions, specifications and requirements.

### Hot Water Requirements

An existing hot water supply may be used as the hot water source for your Glassware Washer. The washer requires the inlet water temperature to be greater than 120°F (49°C) to fully activate powdered detergents and raise the glassware temperature to achieve fast and complete drying. Hotter water will produce better washing results.

Water pressure must be between 20-120 psi (138-827 kPa) at the washer and provide a minimum of 1.25 gallons (4.7 liters) per minute flow rate.

A shut-off valve should be installed in the water supply line plumbed to the washer. The hot water inlet valve on the Glassware Washer is equipped with a female 3/8 NPT fitting. A compression fitting is provided that connects to .375" (9.53mm) OD copper tubing.

### Purified Water Requirements

To use purified water for the rinse cycles, a purified water supply is required. The washer can be connected to an in-house, pressurized, purified water tap; a water purification system; or a non-pressurized purified water container.

The purified water system must be provided with supply piping of sufficient size to permit at least 0.9 gallons (3.4 liters) per minute of flow. Tubing with an internal diameter of at least .375" (9.53 mm) is recommended. The purified water inlet valve on the washer is equipped with a plastic hose barb connection to accommodate .375" (9.53 mm) ID flexible plastic or rubber hose. Use a spring or band hose clamp to secure the hose to the hose barb connection. The hose barb connection may be removed from the valve to expose a male 3/4" - 11-1/2 NH hose coupling. Rigid plastic, tin-lined, or stainless steel tubing and fitting can be connected to the washer fitting.

An optional connection kit (Labconco P/N 4592200) is available for connecting .375" (9.53mm) OD plastic tubing to the purified water inlet valve. The purified water system connections must be airtight to prevent the washer pump from pulling air instead of water.


If the purified water is stored in a carboy, at least 3.4 gallons (13 liters) must be available for each rinse. If all the purified water is not available at the start of a wash cycle, but instead is produced concurrently with the wash cycle, the production rate must permit at least 0.9 gallon (3.4 liters) per minute to be delivered to the washer.

## Electrical Requirements

Undercounter and Freestanding models, are hard-wired directly into a junction box using conduit. A 20 Amp circuit breaker or fuse is required for models rated at 115V (60 Hz) or 230V (50/60 Hz).

High voltage models are designed to operate at 230 VAC, single phase. It will operate safely at nominal 208 VAC, however the heater output will be significantly degraded as the operating voltage drops below 230V. It is highly recommended that if the voltage is low that a boost transformer is used.

If your Glassware Washer is converted to a Mobile model, a dedicated electrical outlet is required. A 20 Amp circuit breaker or fuse is required for models rated at 115V (60 Hz) or 230V (50/60 Hz). 115V models are equipped with a 20 Amp NEMA 5-20P plug. 230V models are equipped with a 15 Amp NEMA 6-15P plug. It may be necessary to remove the plug and install a different plug to match the available receptacle.

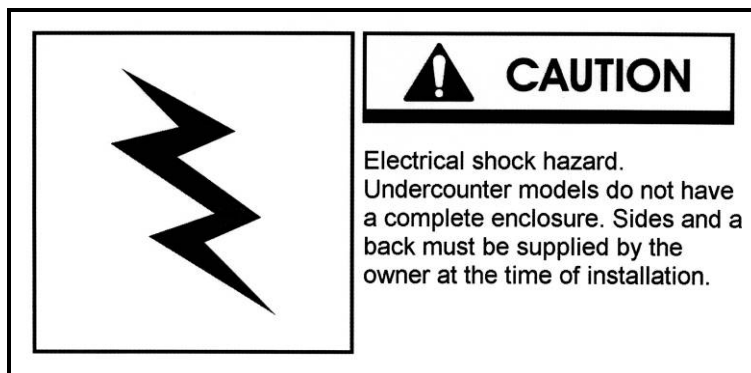



## **Drain Requirements**

The drain hose, .75" (1.9 cm) ID is installed during manufacture. The hose provides a flexible coupling to the building drain piping and can be secured with a spring or band hose clamp. (A band hose clamp is provided with the washer.)

Tubing or pipe, .75" (1.90 cm) ID or larger, should be provided for the building drain. The use of an air gap is strongly recommended to prevent siphoning of wastewater into the washer.

Do not reduce the size of the drain plumbing.

Glassware Washers that have been converted to Mobile models drain into a sink through a hose assembly that attaches to the faucet.

## **Space Requirements**

If your Glassware Washer is an Undercounter model, you must ensure the undercounter opening is the proper size to accommodate the washer. The dimensions for the Undercounter washer are shown in Appendix B.

---

# CHAPTER 3

## GETTING STARTED

Now that the site for your Glassware Washer is properly prepared, you are ready to unpack, inspect, install, and test your washer.

You may need common plumbing and electrical installation tools in addition to a 9/16" wrench, a flat-blade screwdriver, a Phillips screwdriver, and a carpenter level to complete the instructions in this chapter.



**The Glassware Washer weighs over 150 lbs. (68 Kg). The carton allows for lifting with a mechanical lift truck or hand truck. If you must lift the washer manually, use at least two (2) persons and follow safe lifting guidelines.**



**CAUTION**

**Lifting hazard.  
Do not lift or move this  
equipment without  
assistance.**

## Unpacking Your Glassware Washer

Carefully unpack your Glassware Washer and inspect it for damage that may have occurred in transit. If your Glassware Washer is damaged, notify the delivery carrier immediately and retain the entire shipment intact for inspection by the carrier.

The United States Interstate Commerce Commission rules require that claims be filed with the delivery carrier within fifteen (15) days of delivery.

Do not discard the carton or packing material for your Glassware Washer until you have checked all of the components and installed and tested the Glassware Washer.

**NOTE: Do not return goods without the prior authorization of Labconco. Unauthorized returns will not be accepted. If your Glassware Washer was damaged in transit, you must file a claim directly with the freight carrier. Labconco Corporation and its dealers are not responsible for shipping damage.**

Do not remove the washer from its shipping skid until it is ready to be placed into its final location. Move the washer by placing a flat, low dolly under the shipping skid.

## Washer Components

Locate the model of washer you received in the following group of tables. Verify that the components listed are present and undamaged.

| <b>Catalog #</b> | <b>Washer Description</b>   | <b>SteamScrubber Undercounter Models</b> |
|------------------|---|--|
| 4400320          | SteamScrubber Undercounter – 115 V with upper and lower racks             |  |
| 4400321          | SteamScrubber Undercounter – 230 V with upper and lower racks             |  |
| 4400330          | SteamScrubber Undercounter – 115 V with window with upper and lower racks |  |
| 4400331          | SteamScrubber Undercounter – 230 V with window with upper and lower racks |  |

### Plus the Following Loose Parts:

| <b>Part #</b> | <b>Component Description</b>                                   |
|---------------|--|
| 1879801       | Leveling Feet (4)  |
| 4591100       | User Manual  |
| 4462900       | Screws (2)   |
| 1488800       | Clamp (1)  |
| 4522500       | LabSolutions Powder Detergent with MSDS – 6 ounces (175 grams) |
| 1410700       | Water Connector  |
| 1448700       | Coupling & Sleeve (1)  |

| <b>Catalog #</b> | <b>Washer Description</b>   | <b>SteamScrubber Freestanding Models</b> |
|------------------|---|--|
| 4400420          | SteamScrubber Freestanding – 115 V with upper and lower racks             |  |
| 4400421          | SteamScrubber Freestanding – 230 V with upper and lower racks             |  |
| 4400430          | SteamScrubber Freestanding – 115 V with window with upper and lower racks |  |
| 4400431          | SteamScrubber Freestanding – 230 V with window with upper and lower racks |  |

**Plus the Following Loose Parts:**

| <b>Part #</b> | <b>Component Description</b>                                   |
|---------------|--|
| 1879801       | Leveling Feet (4)  |
| 4591100       | User Manual  |
| 4468000       | Hole Plug (2)  |
| 1488800       | Clamp (1)  |
| 4522500       | LabSolutions Powder Detergent with MSDS – 6 ounces (175 grams) |
| 1410700       | Water Connector (1)  |
| 1448700       | Coupling & Sleeve (1)  |

**Catalog # Washer Description**

|         |  |
|---------|--|
| 4420320 | FlaskScrubber Undercounter – 115 V with lower spindle rack             |
| 4420321 | FlaskScrubber Undercounter – 230 V with lower spindle rack             |
| 4420330 | FlaskScrubber Undercounter – 115 V with window with lower spindle rack |
| 4420331 | FlaskScrubber Undercounter – 230 V with window with lower spindle rack |

**FlaskScrubber  
Undercounter  
Models**

**Plus the Following Loose Parts:**

| <b>Part #</b> | <b>Component Description</b>                                   |
|---------------|--|
| 4584600       | Lower Spindle Rack with 36 spindles                            |
| 1879801       | Leveling Feet (4)  |
| 4591100       | User Manual  |
| 4462900       | Screws (2)   |
| 1488800       | Clamp (1)  |
| 4522500       | LabSolutions Powder Detergent with MSDS – 6 ounces (175 grams) |
| 1410700       | Water Connector (1)  |
| 1448700       | Coupling & Sleeve (1)  |
| 4525200       | Spindle – Small (10)   |
| 4414703       | Clip – Spindle (10)  |
| 4425101       | Plug – Spindle (36)  |


**Catalog # Washer Description**

|         |  |
|---------|--|
| 4420420 | FlaskScrubber Freestanding – 115 V with lower spindle rack             |
| 4420421 | FlaskScrubber Freestanding – 230 V with lower spindle rack             |
| 4420430 | FlaskScrubber Freestanding – 115 V with window with lower spindle rack |
| 4420431 | FlaskScrubber Freestanding – 230 V with window with lower spindle rack |

**FlaskScrubber  
Freestanding  
Models**

**Plus the Following Loose Parts:**

| <b>Part #</b> | <b>Component Description</b>                                   |
|---------------|--|
| 4584600       | Lower Spindle Rack with 36 spindles                            |
| 1879801       | Leveling Feet (4)  |
| 4591100       | User Manual  |
| 4468000       | Hole Plug (2)  |
| 1488800       | Clamp (1)  |
| 4522500       | LabSolutions Powder Detergent with MSDS – 6 ounces (175 grams) |
| 1410700       | Water Connector (1)  |
| 1448700       | Coupling & Sleeve (1)  |
| 4525200       | Spindle – Small (10)   |
| 4414703       | Clip – Spindle (10)  |
| 4425101       | Plug – Spindle (36)  |

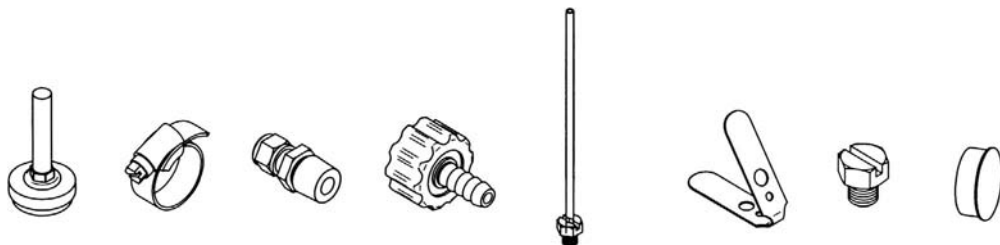

1879801 148880 1410700 1448700 4525200 4414703 4425101 4468000

If you did not receive one or more of the components listed for your washer, or if any of the components are damaged, contact Labconco Corporation immediately for further instructions.

## Removing the Shipping Skid

After you verify the washer components, move your washer to the location where you want to install it. Then, follow the steps listed below to remove the shipping skid from your washer.

### **To remove the shipping skid:**


1. Remove accessories and loose items from inside the washer. Cut tie wraps and remove racks.
2. Gently place the washer on its back, positioning the washer on the shipping carton or other cushioning material.
3. Remove the four screws fastening the skid to the washer.
4. Retain the shipping skid until you test the various features of the washer.

## Washer Installation

After you remove the skid from your washer, you must install leveling feet on the bottom of the washer. If you have an Undercounter or Freestanding model, then you must level the washer and center the washer door. If you have an Undercounter model, you must also secure the washer to the underside of the countertop or worksurface.

## Installing the Leveling Feet


1. Locate the holes on the bottom of the washer where the four skid-fastening screws were previously attached.
2. Thread the four leveling feet into the holes.
3. Carefully move the washer into an upright position. Be careful to distribute the weight evenly on the leveling feet when setting the washer upright.
4. Reinstall racks.

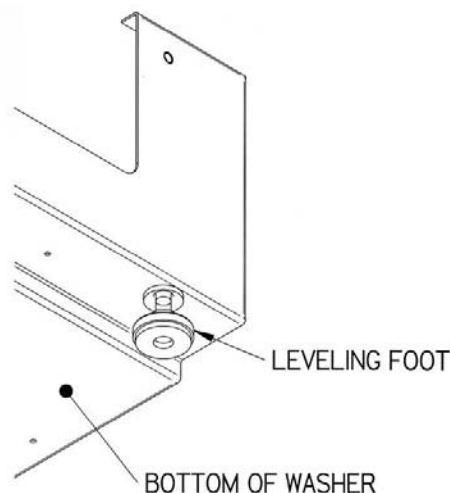


## Leveling the Washer

If you have a Freestanding model of the SteamScrubber or FlaskScrubber Glassware Washer, after you have installed the leveling feet, level the washer by adjusting the leveling feet.


1. Position the washer in its final installed location.
2. Place a carpenter level on the washer in the positions shown.

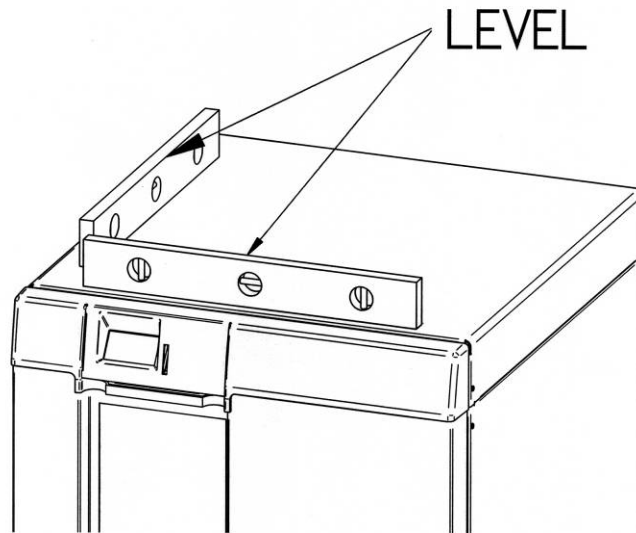

3. Turn the four leveling feet, as needed, to level the washer.

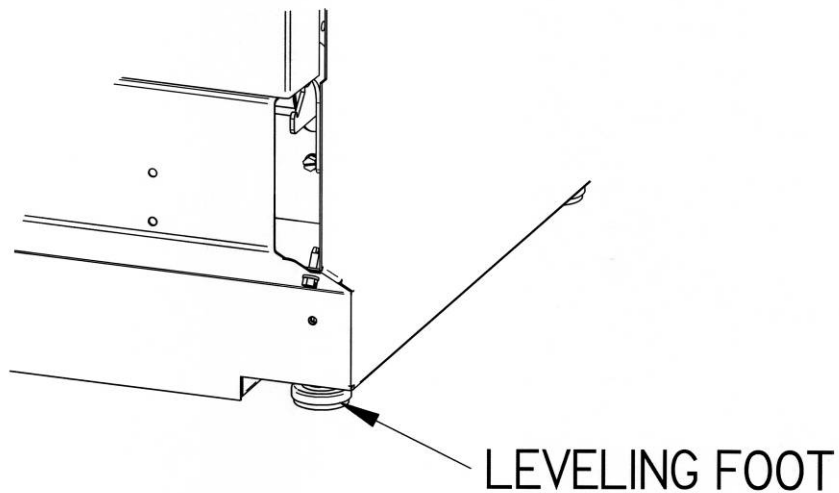

If you have an Undercounter model, level it side-to-side by placing a short level on the inside top edge of the door. See the figure below for the placement of the level on the washer. Check the level front to back by placing a level on the rail that the lower rack slides on.

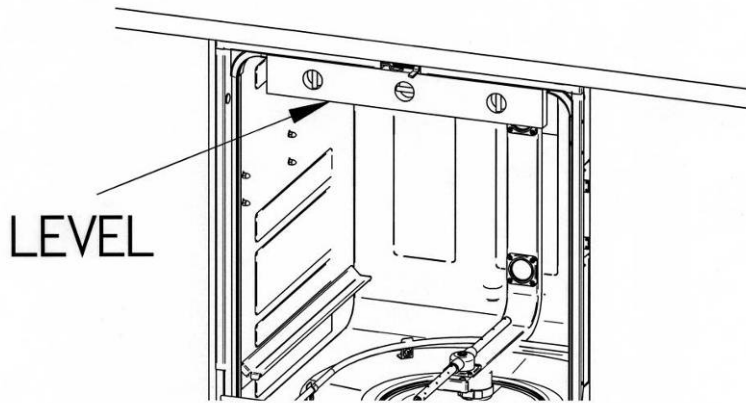


### Centering the Washer Door

When the washer is level, check the operation of the washer door for alignment with the washer tank. If the door is not centered in the opening, you must center the door by adjusting the corresponding leveling foot. (For example, if the door hits the right side of the tank, raise the right front corner of the washer. If the door hits the left side of the tank, raise the left front corner of the washer.)

### Securing the Washer to the Countertop

If you have an Undercounter Glassware Washer, two screws have been provided to attach the front of the washer to the underside of the countertop or work surface. After the washer is leveled in its final position and the washer door is properly centered, install the two screws as shown in the figure below to stabilize the unit and prevent it from moving.

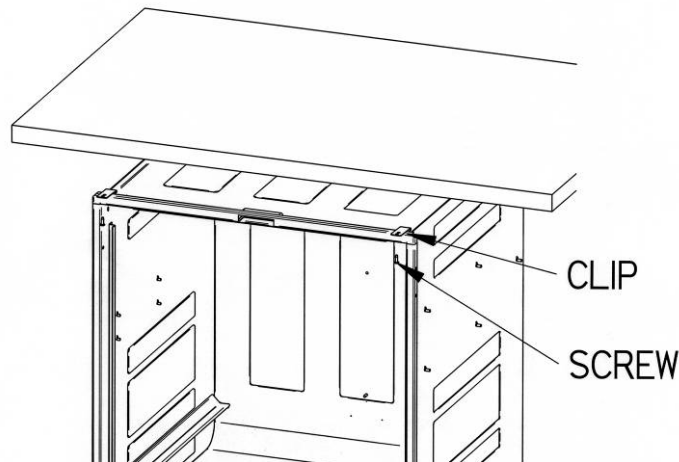

## Connecting the Water Services

A hot water supply must be attached to your Glassware Washer. In addition, a purified water source must be connected to your washer if pure water rinses will be required.


## Connecting the Hot Water Supply

To prevent valve clogging, thoroughly flush the hot water lines for your hot water supply source prior to connecting the washer to the water lines. The water supply valve to which the hot water supply is connected is located on the back of the washer base. The back panel of FreeStanding models must be removed to access the valve.

You must provide the plumbing fittings and components needed to connect the hot water supply source. The washer hot water supply valve is equipped with a compression fitting for a .375" (9.53 mm) OD tube.

### To connect the hot water supply:

1. Be certain the hot water supply piping has been thoroughly flushed prior to connecting it to the washer. Debris in the piping can clog the washer valve.
2. Make the necessary plumbing connections to properly connect the hot water supply to the washer.
3. It may be necessary to place a loop in the supply tube to allow the washer to be slid in or out of its operating position. Be careful not to kink the water supply tube when the washer is placed into its final location. If kinking occurs, it may be necessary to replace the straight fitting on the valve with an elbow.

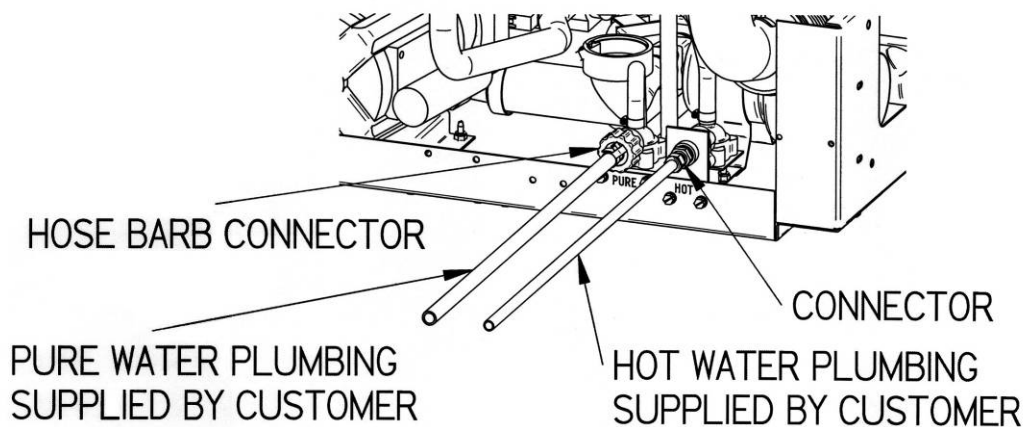

Washer Back View



## Connecting the Purified Water Supply

To prevent valve and pump clogging, flush all of the water lines for your purified water supply source prior to connecting the washer to the water lines. The valve to which the purified water supply must be connected is located on the back of the washer base. The back panel of Freestanding models must be removed to access the valve. Refer to the previous figure.

The purified water inlet valve is equipped with a plastic hose barb connection to accommodate .375" (9.53 mm) ID flexible plastic or rubber hose. The hose connection may be removed from the valve to expose a male 3/4" – 1 1/2 NH hose coupling. Rigid plastic, tin-lined, or stainless steel tubing and fittings may be connected to the washer fitting.

At least 0.9 gallon (3.4 liters) per minute of purified water flow is required. At least 3.4 gallons (12.9 liters) of purified water must be available for each pure rinse. A .375" (9.5 mm) ID pipe is the minimum recommended.

If you do not use purified water during the rinse cycles, you may disable the purified water pump. Refer to Controller Set-Up in this section.


If the washer has been converted to a Mobile model and is moved to various locations, and if purified water is used, provisions must be made to use a flexible hose from a source of house-purified water. If a carboy is used, place the carboy on a carboy cart and move it with the washer. Be careful not to let the hose kink or become damaged.

### **To connect the purified water supply:**


1. Make certain the purified water supply piping has been thoroughly flushed prior to connecting it to the washer.
2. Make the necessary plumbing connections to properly connect the purified water supply to the washer.

**NOTE:** If the purified water source is a carboy, complete steps 3 and 4.

3. Connect the purified water source through either the top or the bottom of the carboy, as illustrated in the following figures. If the purified water supply feeds directly from the carboy top, make certain the hose is submerged under the water line. Attach a weight to the end of the tubing to keep it from floating on top of the water and possibly allowing air into the system.
4. Make certain the top of the carboy is vented to permit proper flow from the carboy to the washer. Make certain that all of the connections on the purified water system are air tight. The pump must not pull air instead of water. Make certain the supply hose is not crimped.

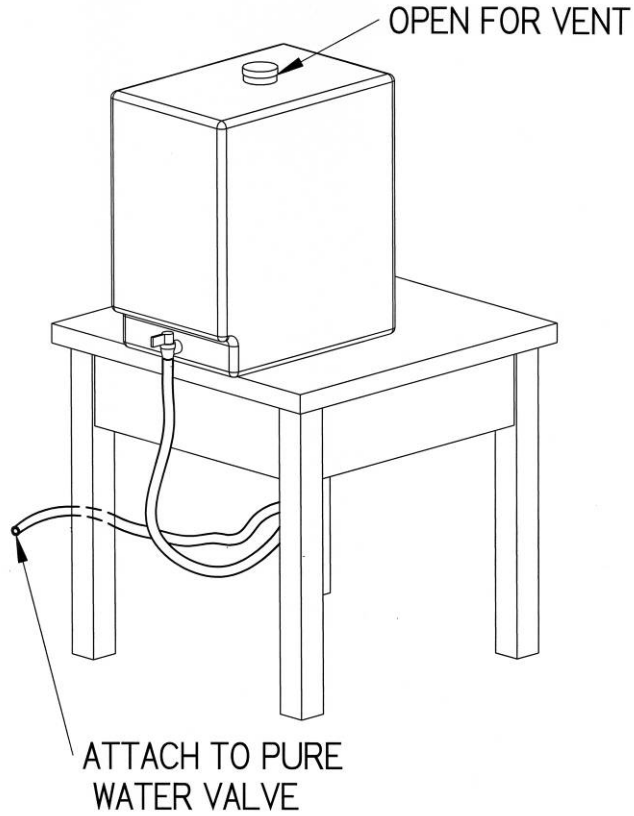

**Example of a Bottom Feed Connection**

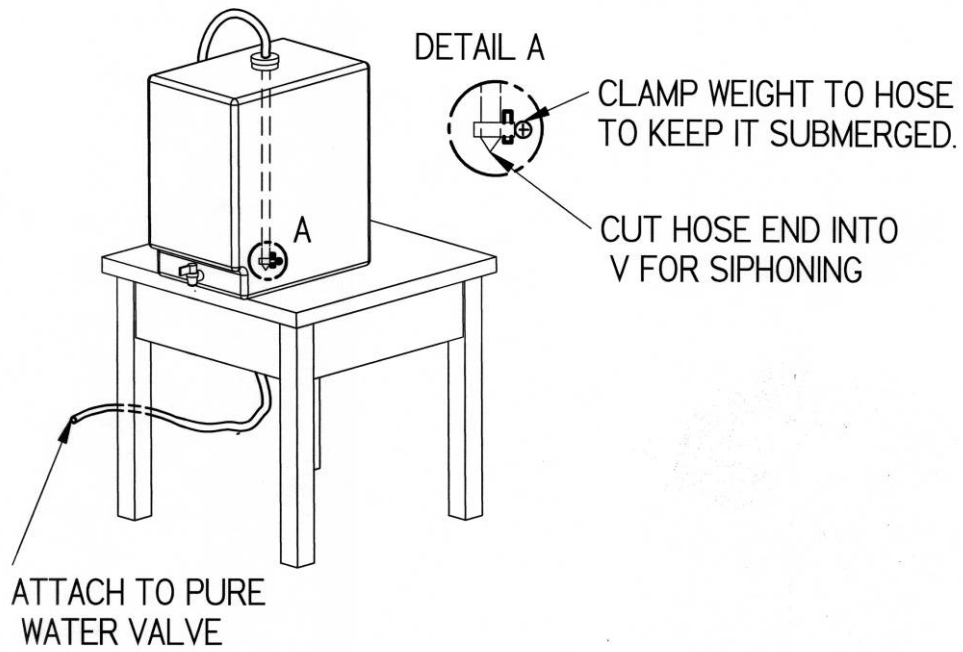

**Example of a Top Feed Connection**

## Connecting the Drain Hose

Before continuing with this section, be certain that your drain meets the requirements detailed in *Drain Requirements* in *Chapter 2: Prerequisites*.

Be certain you conform to all local plumbing codes when you route and connect the drain.

The drain hose is installed at the factory. The hose provides a flexible coupling to the drain piping and can be secured with a spring or band hose clamp. (A band hose clamp is provided with your washer.)

Tubing or pipe .75" (1.9 cm) ID or larger should be used for the drain.


Do not route the drain to an elevation more than 30" (76 cm) above the pump discharge. Do not reduce the size of the drain plumbing.

Do not remove the hose from the clamp on the rear of the glassware washer.

The use of an air gap is strongly recommended for all installations to prevent the siphoning of wastewater into the washer.


The drain piping can be routed to either a sink or a floor drain. If you route the drain piping to a sink drain without using an air gap, the highest point of the drain pipe must be above the highest water level in the sink, as shown below.

If an air gap is not used, the drain line must be elevated above the water level of the sink.

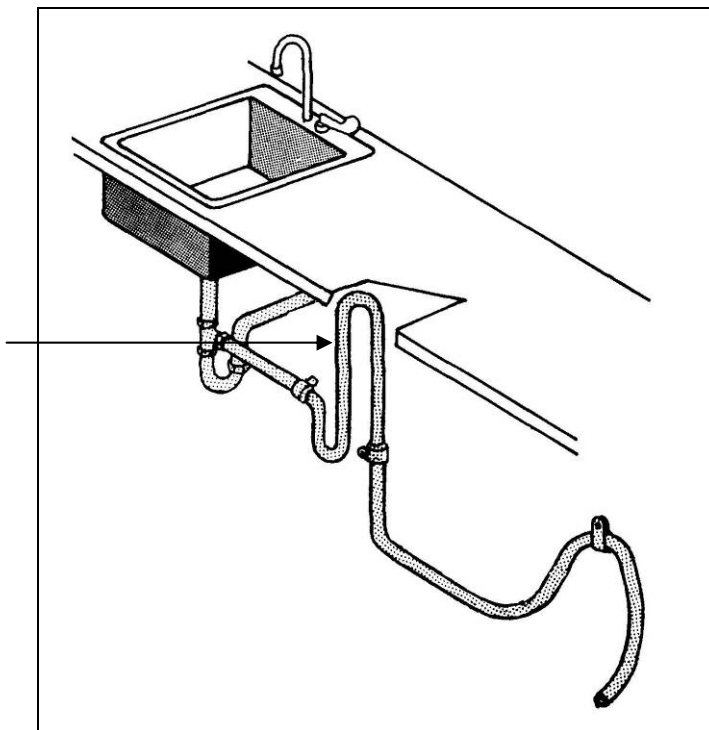


### Chapter 3: Getting Started

If you route the drain piping to a sink drain and use an air gap, refer to the figure below for the placement of the air gap.

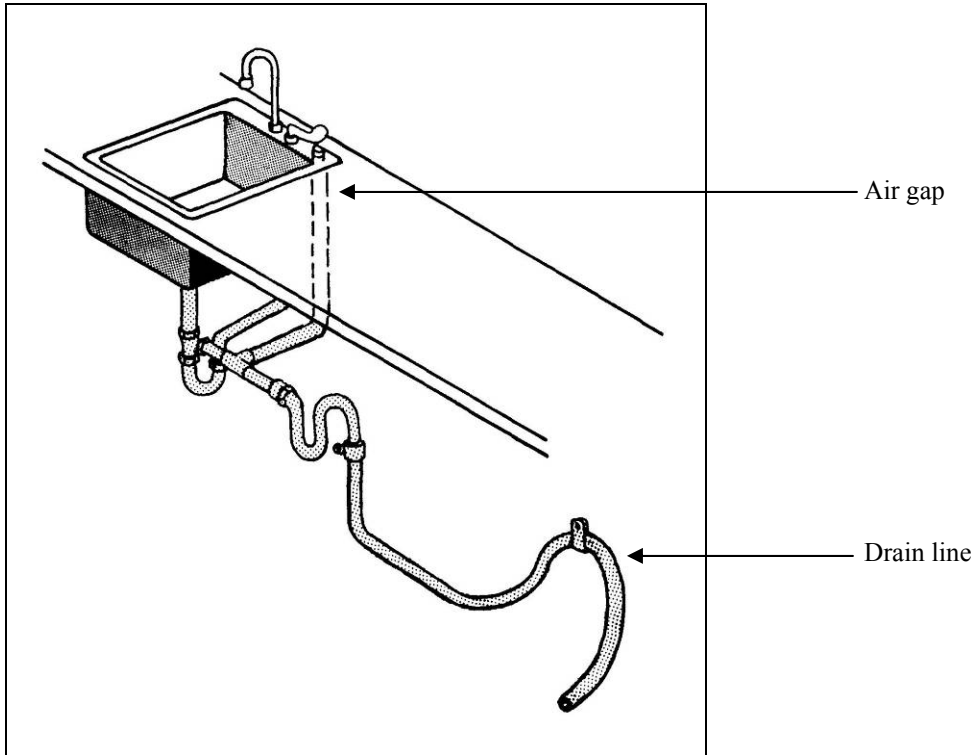

If you route the drain piping to a floor drain, refer to the figure below for placement of the drain tubing and trap.

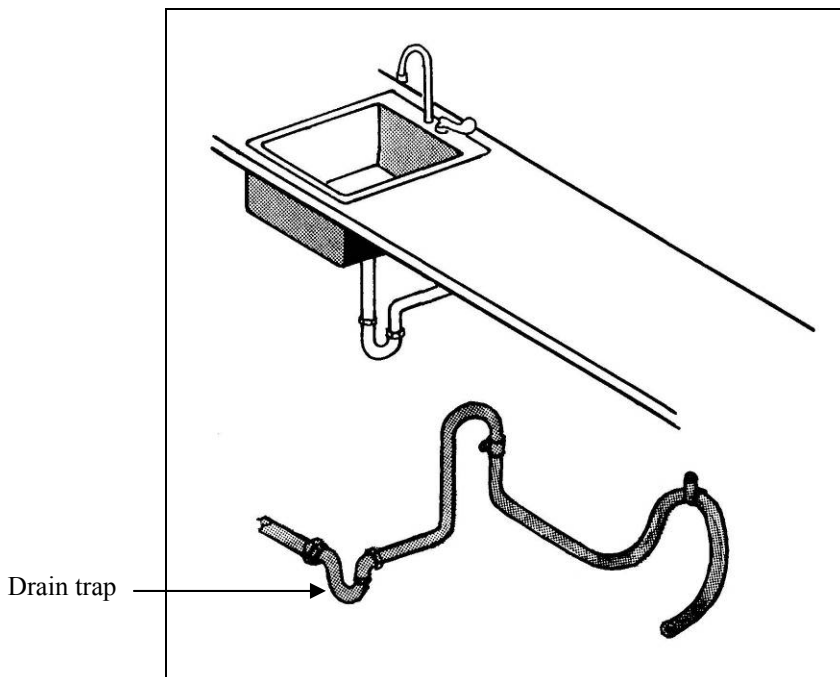

## Connecting the Power Supply

Before continuing with this section, be certain your power supply meets the requirements specified on the washer data plate and in *Electrical Requirements* in *Chapter 2: Prerequisites*. Undercounter and Freestanding Washers should be hard-wired directly into a junction box. Refer to the figure below and follow the steps below to connect your washer to the appropriate power source.


Be certain that you check and conform to all local electrical codes. Do not disturb any internal wiring in the washer.

If you are connecting to certain legs of a 3-phase power system to obtain single-phase power, be sure to consult a professional to insure codes/standards are followed. These professionals can determine the appropriate connections for either a Wye or Delta connected power source.

The supply circuit must consist of number 12 AWG 3-wire cable or heavier. Washers must be connected to a dedicated service with a 20 amp circuit breaker or fuse. Wire must comply with all local electrical codes. Type THW or THWN wire may be used if it complies.


### **To connect the electrical supply:**

1. Remove the lower panel on the washer front by removing the two screws located at the bottom of the panel.

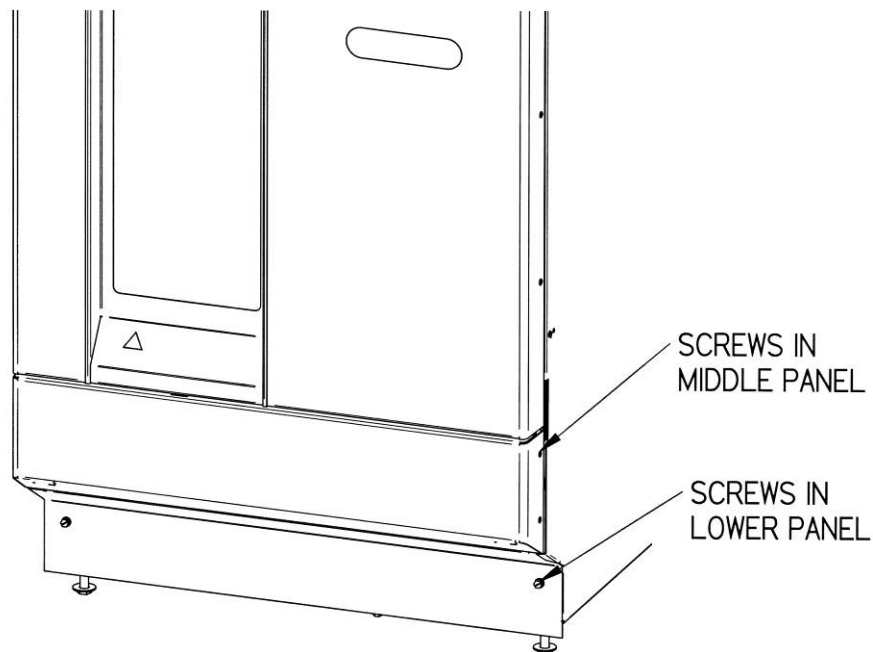

2. Remove 2 screws on each side of the middle panel.
3. Pull the middle panel forward and remove it.


- Remove the cover over the electrical junction box by removing the two screws.

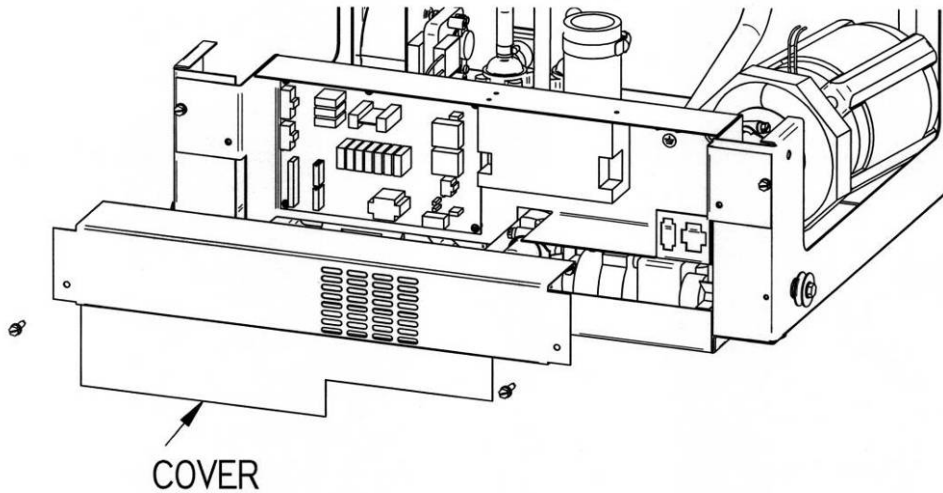

- Make sure electrical power is turned off before making any connections.
- On 115V models, connect the customer-supplied power wires to the black and white wires, using wire connectors. On 230V models, attach customer-supplied power wires to the brown and blue wires, using wire connectors. Customer-supplied wires must be clamped to the hole where they pass through the rear of the junction box, if prescribed by local codes. If it is necessary to use a bushing with an integral clamp at the inlet to the junction box, remove the plastic bushing supplied with the washer before installing the new customer supplied bushing.
- Connect the customer-supplied ground wire to the green screw.
- Replace the cover on the electrical junction box.
- Replace the lower panel and door panel assembly on the washer.

## Controller Set-Up

The washer control system has been set at the factory to agree with the configuration of the washer as it was shipped. The control may be customized or modified if certain optional accessories are added.

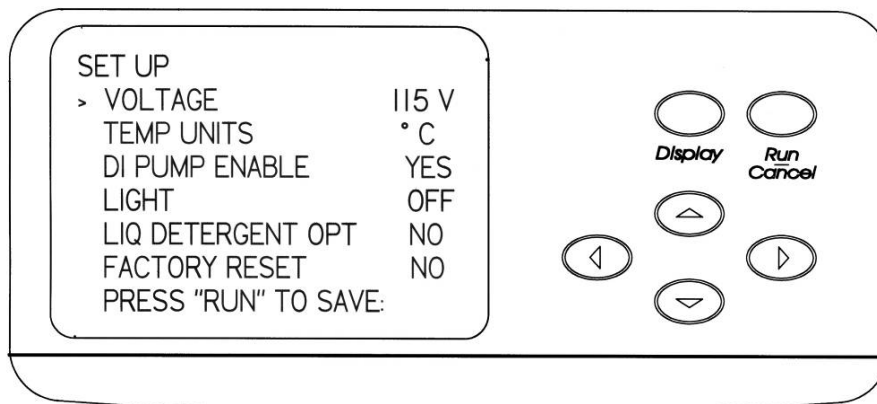

**To set up the washer control:**


1. Locate the control panel on the front of the washer.
2. Press and hold the ▼ button and close the handle of the washer by sliding the door latch knob to the right. Hold for 5 seconds.
3. The display will show the SET UP screen.
4. The > on the left side will highlight VOLTAGE.
5. Press the ◀ or ▶ buttons to select the correct operating voltage of the washer.  
CAUTION: The voltage selected (115V or 230V) does not change the acceptable AC mains voltage that can be connected to the washer. It only selects the programs that are suitable for the identified voltage.
6. Press the ▼ button and TEMP UNITS will highlight.
7. Press ◀ or ▶ buttons to select either °F or °C which will be used on the display to set and monitor temperature.
8. Press the ▼ button and DI PUMP ENABLE will highlight.
9. Press ◀ or ▶ buttons to YES or NO. This allows the pure water valve and pump to be disabled if pure water will not be used for rinse cycles and prevents inadvertently programming a cycle to have pure rinses.
10. Press the ▼ button and LIGHT will be highlighted on the display.
11. Press ◀ or ▶ to cause the light inside the tank on window models to be turned on or off while the door is latched.
12. Press the ▼ button and LIQ DETERGENT OPT will be highlighted on the display.
13. Press ◀ or ▶ to activate the control for the liquid detergent dispense option if this accessory has been installed.
14. Press the ▼ button and FACTORY RESET will be highlighted on the display.
15. Press the ◀ or ▶ to select YES to erase all program set points entered by the user and to restore all set points that were active when the washer was shipped.
16. Press RUN to save all settings.

**Testing Your Washer**

Now that you have made the required water and electrical connections to your washer, a series of diagnostic steps are provided in this section to test the operation of the washer.


Before turning on the washer, make certain the water is turned on and no packing material remains inside the washer.

If the washer is interrupted in the middle of the diagnostic steps, wait 4 to 5 seconds before opening the washer door to prevent hot water from splashing out. The hot water and steam from the washer may cause scalding.

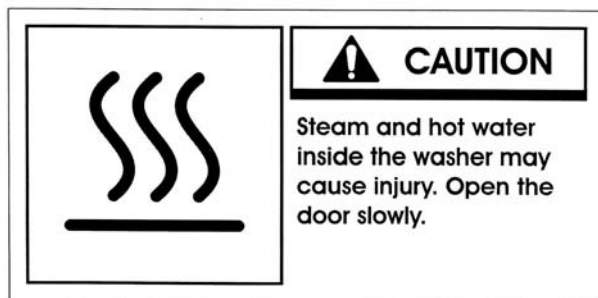

## Running the Diagnostics

To ensure your washer is operating properly, perform the diagnostics procedure detailed below, which will take approximately 10 minutes to complete. Refer to the figure below to familiarize yourself with the control panel before you begin the procedure.

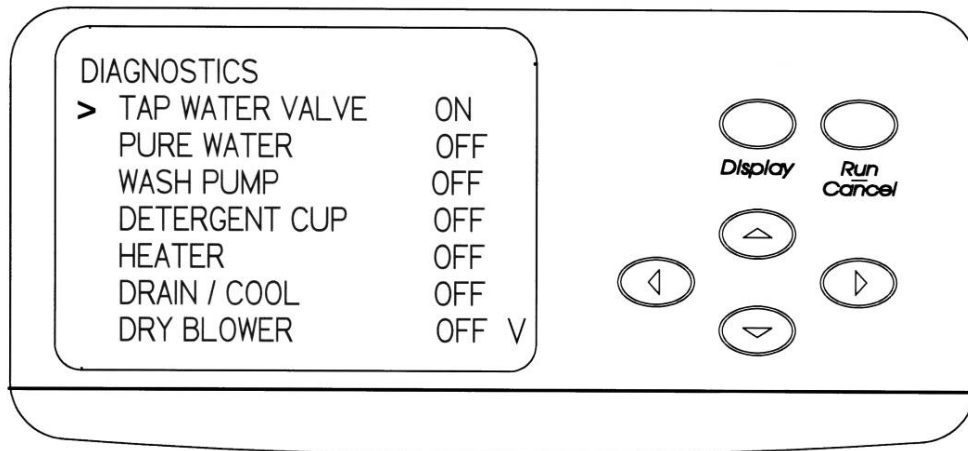

### To run the diagnostics test:

1. Press the ▲ button on the control panel and close the door and slide the door latch to the right. Hold for 5 seconds.
2. The display will show the Diagnostics screen and the indicator on the left side will point to TAP WATER VALVE.
3. Press either the ► or ◀ button to turn the tap water valve ON. Tap water should enter the washer. Allow water to enter for approximately 1 minute. Press either the ► or ◀ button to turn the water off.
4. Press the ▼ button and the arrow on the display will point to PURE WATER.
5. Press either the ► or ◀ button turn the pure water pump ON and open the pure water valve. Allow pure water to enter the washer. After about 3 minutes the fill switch should cause the water flow to turn off.
6. Press the ▼ button and the arrow on the display will point to WASH PUMP.
7. Press the ► or ◀ button and the wash pump should energize.
8. Press the ▼ button and the arrow on the display will point to DETERGENT CUP.
9. Press either the ► or ◀ button to energize the detergent cup. Within about 1 minute the detergent cup door should be heard to open.
10. Press the ▼ button and the arrow on the display will point to HEATER.
11. Press either the ► or ◀ button and the heater will be energized.
12. Press the ▼ button and the arrow on the display will point to DRAIN / COOL.
13. Press either the ► or ◀ button and the drain will be energized. Allow all water to be drained from the washer. The sound of the pump will change when all water is drained. This will take about 1-1/2 to 2 minutes. If the Cool Water Drain Kit accessory has been installed, cool water will mix with hot water draining from the washer.



14. Press the ▼ button and the arrow on the display will point to DRY BLOWER.
15. Press either the ► or ◀ button and the dry cycle blower will be energized.
16. The ▼ arrow in the lower right side of the display indicates that more selections may be made. Press the ▼ button to select the next function to test. The arrow on the display will point to LIQ DETERG PUMP.
17. Press the ► or ◀ button and if the liquid detergent accessory has been installed and if it was activated in the Set Up procedure, the detergent pump will operate.
18. Press the ▼ button and the arrow on the display will point to LIGHT.
19. Press either the ► or ◀ button and if the washer is equipped with a window in the door, the light inside the washer will illuminate.
20. Press the ▼ button and the arrow on the display will point to COOL FAN.
21. Press either the ► or ◀ button and the small fan inside the kick panel will operate.
22. This completes the check of most of the washer functions. Unlatch the door to exit from the Diagnostics mode.
23. To verify the proper operation of the overfill switch, repeat steps 1 & 2 above. Press either ► or ◀ button and allow tap water to enter the tank until the water flow turns off.
24. Press and hold the RUN button. Water will enter the washer and should turn off before it overflows.
25. Press the ▼ arrow until DRAIN/COOL is highlighted.
26. Press ► or ◀ button to drain water from the washer.
27. Unlatch the door to exit from the Diagnostics mode.

---

# CHAPTER 4

## USING YOUR WASHER

After your washer has passed the basic diagnostics test detailed in *Chapter 3: Getting Started*, you are ready to begin using your washer.

### Arranging the Washer Racks

The FlaskScrubber and SteamScrubber are equipped with different rack configurations to accommodate the various types of laboratory glassware. In addition to the standard configuration of the Glassware Washers, many optional racks and inserts are available to customize your washer to best suit your needs. For a complete list of available rack accessories, refer to *Appendix A: Glassware Washer Accessories*.


If you purchased a SteamScrubber, proceed to the following section, *The SteamScrubber Racks*. If you purchased a FlaskScrubber, proceed to the section, *The FlaskScrubber Racks*, later in this chapter. A FlaskScrubber can be converted to a SteamScrubber by replacing the lower rack and adding an upper rack and slides. Drying will be less efficient than a SteamScrubber. A spindle rack may be used in a SteamScrubber. The interior of the glassware may not dry well.

### The SteamScrubber Racks

The SteamScrubber is equipped with a top rack and a bottom rack. Both racks are designed to accommodate inserts that hold flasks, beakers, culture tubes, graduated cylinders, Erlenmeyer flasks, petri dishes, watch glasses, BOD bottles, and a variety of other laboratory glassware. These inserts are sold separately and are detailed in *Appendix A*.

## Removing the Upper Rack

The upper rack may be easily removed from the Glassware Washer to facilitate loading the rack outside the washer or to allow very tall glassware to be loaded and washed on the lower rack.

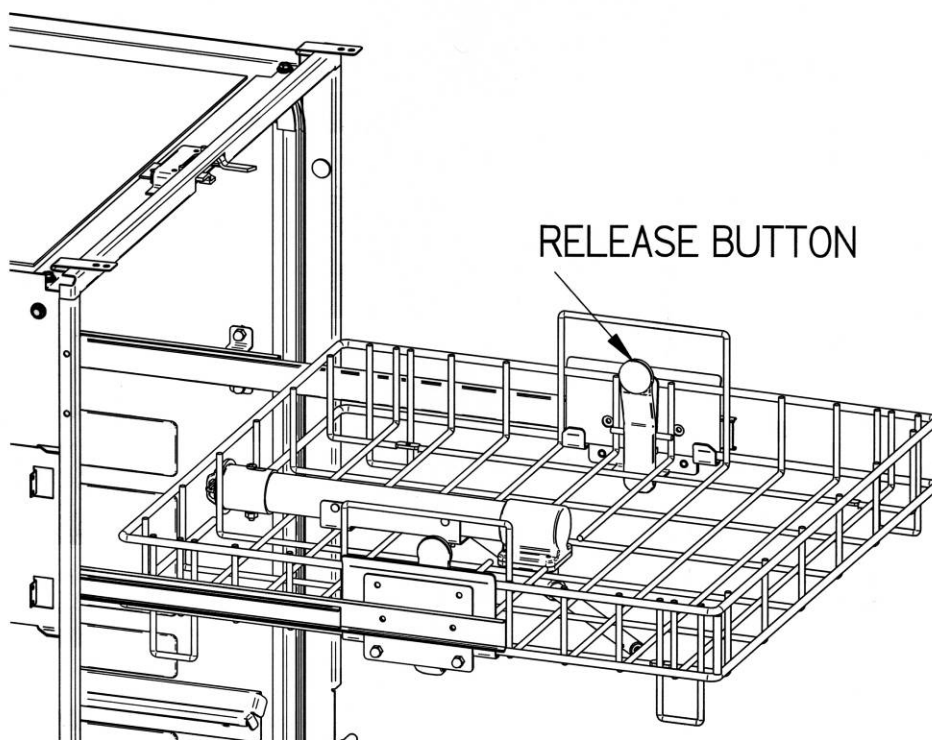

### To remove the upper rack:

1. Slide the upper rack all the way out of the washer.
2. Grasp each side of the rack placing a thumb on the release buttons.
3. Press each release button and lift the rack straight upward.

### To install the upper rack:

1. The upper rack can be replaced into the washer by first sliding each slide rail all the way out of the washer and then lowering the rack onto the rails.
2. Press down and the latch will automatically engage the rail.

## The FlaskScrubber Spindle Rack


The FlaskScrubber Glassware Washer is equipped with a lower spindle rack and glassware holders. The FlaskScrubber spindle rack is designed for narrow-necked glassware such as volumetric flasks, Erlenmeyer flasks, graduated cylinders, Kjeldahl flasks, test tubes, and digestion tubes.

Two sizes of spindles are provided with the washer: thirty-six .25" (6.35 mm) OD and ten .125" (3.18mm ) OD spindles. The optional Forty-Eight Pin Insert (Labconco P/N 4591601) may be positioned directly over spindles or spindles can be removed with a screw driver and replaced with the plugs included with the washer.

Optional top racks are available for the FlaskScrubber to expand the washer's capacity to accommodate many types of glassware. An Upper Spindle Rack (Labconco P/N 4595600) is available with 30 spindles.

### Loading Glassware into the FlaskScrubber

Glassware Holders are supplied with each FlaskScrubber. These holders enable you to position glassware onto the FlaskScrubber spindles.

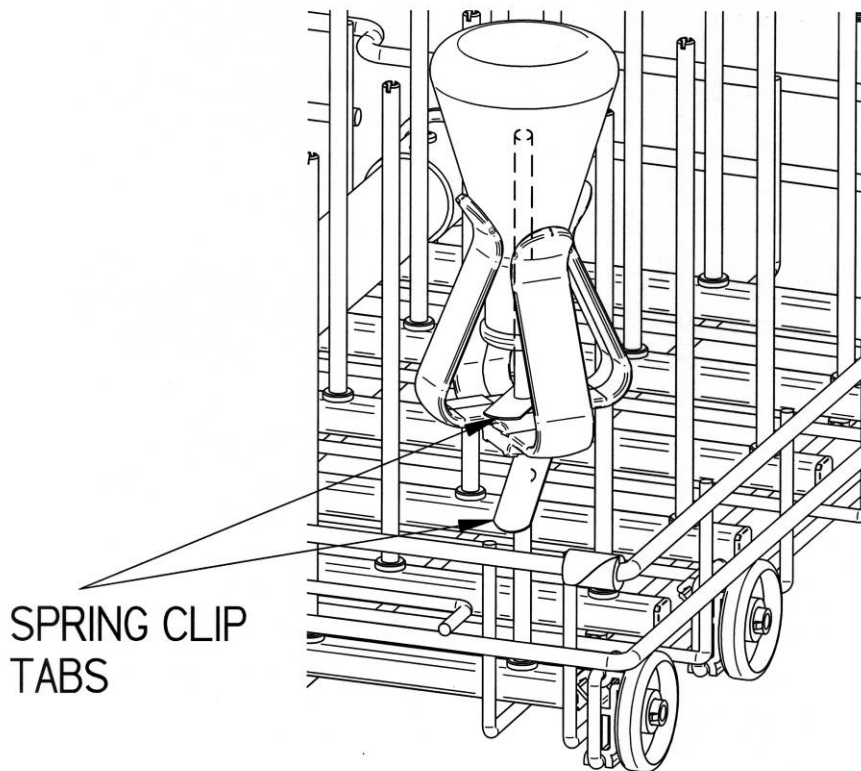


#### **To use a Glassware Holder:**

1. Press the glassware into the holder; the clamps should spring out and grip the glassware.  
**Note:** It is normal for the clamps to deform if large glassware is inserted into the Holder and the washer is operated at high temperature.
2. The holder may be raised or lowered by squeezing the spring clip on the bottom of the holder and then sliding the holder up or down.
3. For best results, do not allow the glassware to rest directly on top of the spindle.

Glassware should be loaded so that it does not touch during a wash cycle. The washer has been designed for quiet operation. To prevent noise and glass breakage, be certain that glassware is retained.

## The Detergent Dispenser

Your Glassware Washer is equipped with a detergent dispenser that contains a main wash cavity and a neutralizing solution port, as shown below. In addition, the washer door has a depression where detergent may be added for use during the first wash of the cycle.

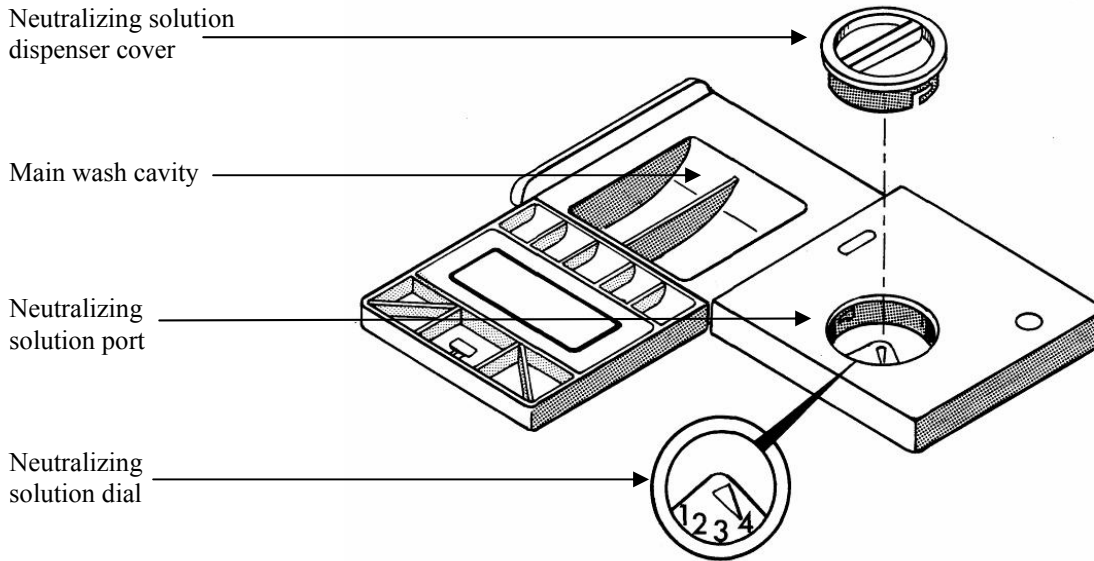


The detergent dispenser opens during the second wash cycle, except when the Plastic cycle is selected, which has only one wash cycle. If your glassware is heavily soiled, you may want to add powder detergent to the depression in the washer door. Do not use liquid detergent in the pre-wash cavity in the washer door. The initial drain and fill cycle flushes the liquid detergent out of the washer prior to the first wash. Refer to the table below to determine which cavities to fill, based on the selected wash choice. You may use either the pre-wash cavity stamped in the washer door or the one on the lid of the detergent cup.

| Wash Cycle       | Fill Pre-wash Cavity | Fill Main Cavity |
|------------------|----------------------|------------------|
| Rinse Only       | NO                   | NO               |
| Plastic          | NO                   | YES              |
| All other cycles | OPTIONAL             | YES              |



**Do not use combustible solvents in the Glassware Washer. If soiled glassware contains flammable compounds, rinse the glassware thoroughly prior to placing the glassware into the washer. Heaters in the washer are very hot and could ignite flammable solvents.**

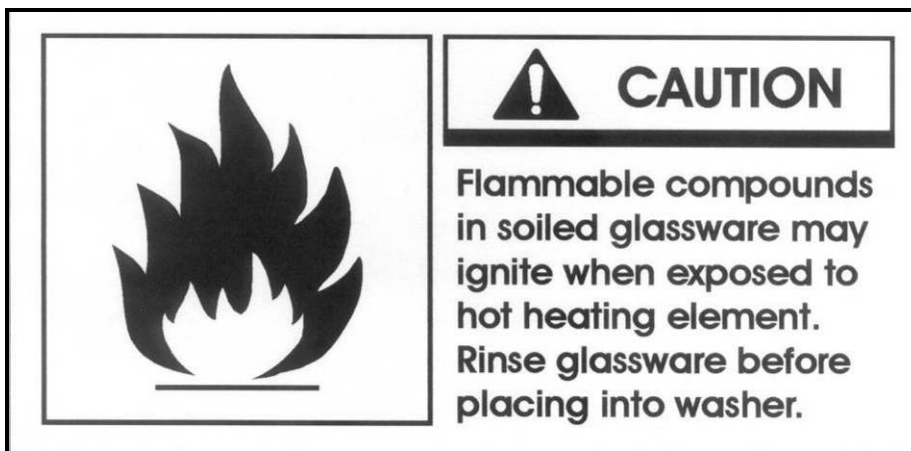

Non-foaming powder or liquid detergent formulated for labware washers, in conjunction with adequate hot water, is essential for clean glassware. Labconco LabSolutions Detergents are recommended. See *Appendix A* for ordering information. Use detergents sparingly. Too much detergent, particularly with softened water may cause etching of the glassware. This appears as a permanent cloudy film on the glassware.

Fill the main detergent cavity three-fourths full if the water hardness grains per gallon (GPG) for your water measures 0-4 GPG (soft water). Fill the detergent cavity completely if your water measures 4 GPG or higher (medium to hard water). If you use the pre-wash cavity stamped in the washer door, fill it completely.

**Note: To open the detergent dispenser, gently slide the white tab sideways. The tab may break if it is pulled upward or perpendicular to the door.**

Store detergent in a cool, dry place. Immediately replace the detergent container lid after dispensing. If lumps develop due to exposure to moisture, discard the detergent according to local regulations.

## Filling the Neutralizing Acid Dispenser


A mild or weak acid may be dispensed from the neutralizing solution port to neutralize the detergent alkalinity. Labconco LabSolutions Neutralizing Acid Rinse is recommended. See *Appendix A* for ordering information. The neutralizing acid rinse is dispensed during the first rinse. To avoid dispensing neutralizing acid during a purified water rinse, the total rinses should be greater than the number of pure rinses. The dispenser holds approximately 170 ml and should be checked and replenished as needed. To fill the neutralizing acid dispenser, refer to the figure above and follow the instructions below.

**To fill the neutralizing acid dispenser:**

1. Twist and pull up on the cap to remove the round cover on the neutralizing solution port.
2. Fill the cavity with LabSolutions Neutralizing Acid Rinse (Labconco P/N 4522200) 1M citric acid or 10% phosphoric acid until the level of the liquid rises and is visible in the port.
3. Set the dial to read 4 ml.
4. Replace the cover.

**Selecting an Operating Cycle**

The microprocessor control on your Glassware Washer allows you to select a factory programmed wash program, to modify the factory preset wash programs or to customize two additional programs. The control panel is shown below.

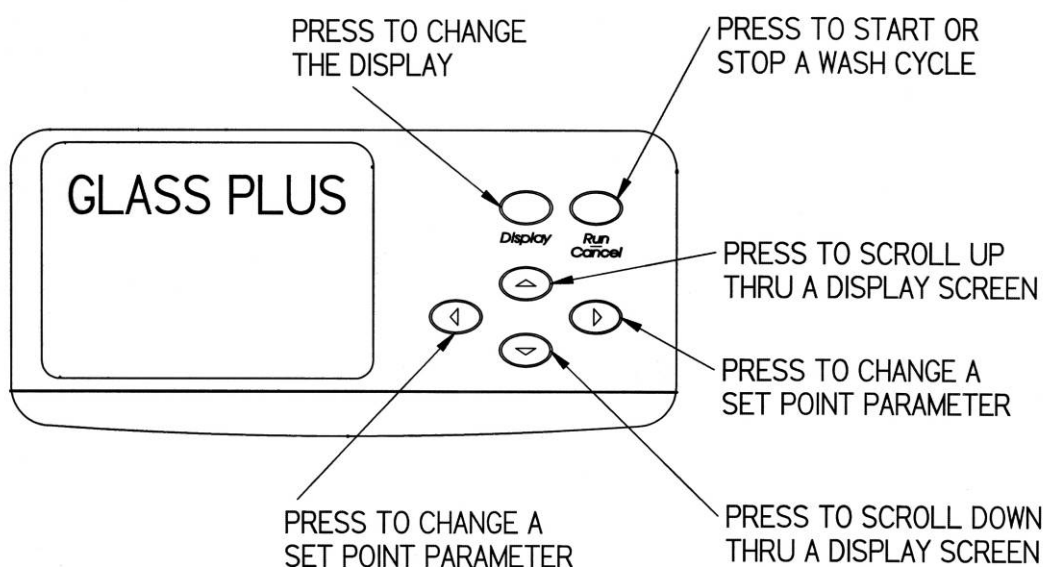**Programming and Starting a Wash Cycle**

Refer to *Appendix C: Glassware Washer Specifications* for complete details on the various factory-set programs and the options available for customizing programs. After you have decided which washer program to use, follow the instructions below to start the washer.


Washers have the capability to wash glassware at high temperatures. Use care when selecting high-heat programs if plasticware is being washed.

Any of the standard factory wash programs may be customized to suit specific wash requirements. All changes made to the programs will be stored in the washer microprocessor's memory until it is changed again. If a FACTORY RESET is performed, all inputs entered by the user will be erased and the programs shipped with the washer will be reinstated. It is possible to customize the following program parameters:

- Wash 2 temperature
- Wash 2 time
- Total number of rinses
- Number of purified water rinses
- Final rinse temperature
- Enable or disable the steam function
- Dry temperature
- Dry time


### **To program and start a wash cycle:**

1. Close the washer door and slide the latch to the right. This energizes the electronic control and the display will show the last program that was run.
2. Press RUN to select this program.
3. If a different program is desired, press DISPLAY. The available programs will be shown on the display.

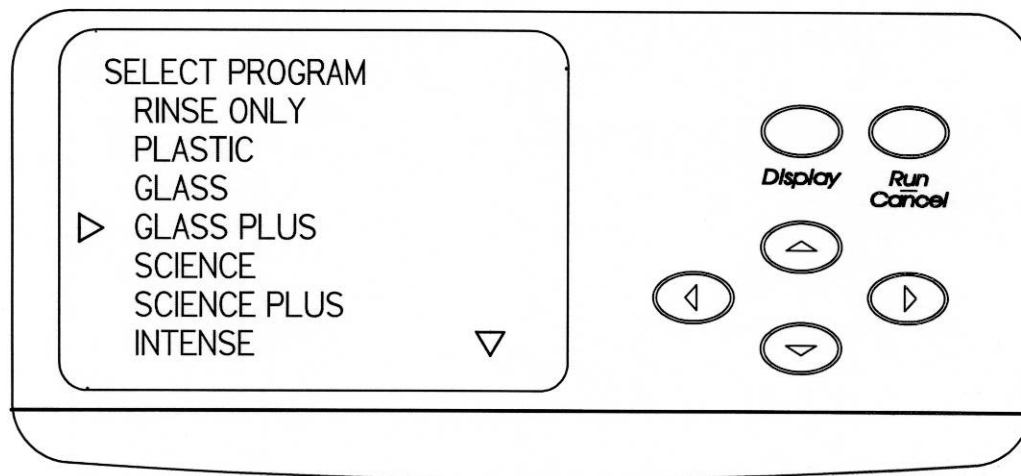

4. Press the ▲ or ▼ buttons until the desired program is indicated by the arrow on the left side. The arrow on the lower right side means that there are additional program selections that will be shown on the display if the ▼ button is pressed. If an arrow is visible in the upper right corner of the display, additional program selections will be displayed when the ▲ button is pressed.
5. Press RUN if no modifications to the program are necessary.
6. If modifications to the program are necessary or you want to see the parameters that are entered into the program, press DISPLAY and the elements in the program that may be altered will be displayed.




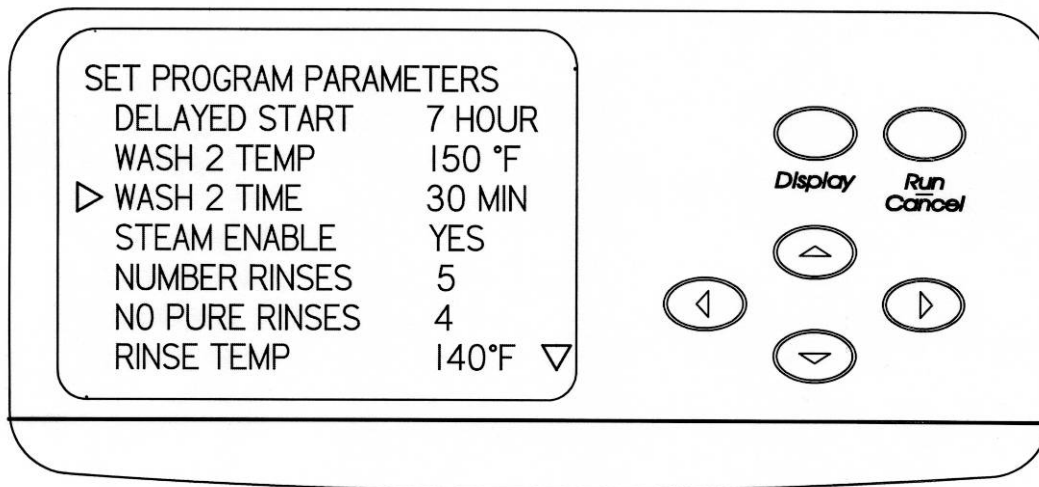

7. Press the ▲ and ▼ buttons to select the parameter that must be altered as indicated by the arrow on the left side. Press ◀ and ▶ buttons to select the appropriate set point. The arrow on the lower right side means that there are additional parameters selections that will be shown on the display if the ▼ button is pressed. If an arrow is visible in the upper right corner of the display, additional parameter selections will be displayed when the ▲ button is pressed.
8. Press RUN when all parameters are satisfactory.

## Interrupting a Wash Cycle

To temporarily pause a wash cycle, simply unlatch the washer door. After the door is unlatched, wait 4 to 5 seconds before opening the door to prevent water from splashing out of the washer.

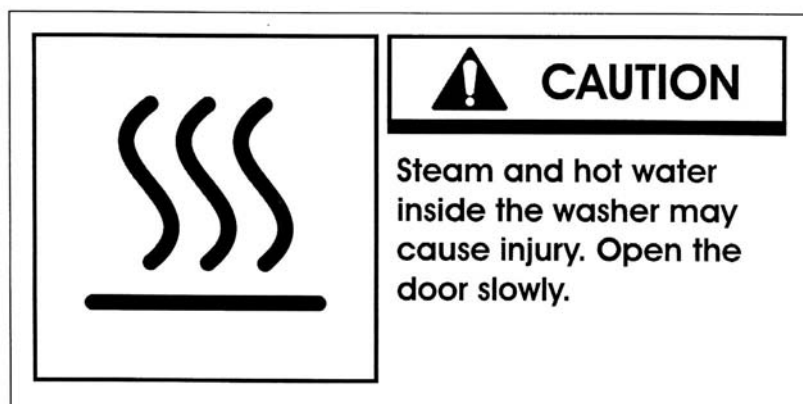

To resume the wash cycle, close and re-latch the washer door. The washer will pause to equalize temperatures and then the cycle will resume at the point it was interrupted.

The wash cycle can be terminated at any time by pressing the Start/Cancel button.

## Alarms

Throughout a wash cycle there are several abnormal events that could occur. The washer is programmed to monitor these occurrences and provide an alarm to alert the operator. Information is shown on the display to identify the alarm situation. The following messages may be displayed. Certain alarm situations are signaled with an audible sound which self cancels after 1 minute.

### HOT GLASS

This displays at the end of a cycle if the temperature inside the washer is over 122°F (50°C). The alarm cancels when the temperature cools.

### PLEASE WAIT

This displays if the door is opened and then closed during a wash or rinse segment of a cycle. The display will also show the time remaining until the operation of the washer will resume where it was interrupted.

### WATER LOW

This displays if a water fill is inadequate. The washer will cancel the cycle and drain water that is in the washer. The alarm is canceled by opening the door. See *Chapter 7: Troubleshooting – Washer will not fill.*

### WATER HIGH

This displays if the washer overfills with water. Water fill valves will be closed and the washer will cancel the cycle and drain water that is in the washer. The alarm is canceled by opening the door. See *Chapter 7: Troubleshooting – Washer overfills.*

### OVERFLOW

This displays if water is detected in the base of the washer. Water fill valves will be closed and the washer will cancel the cycle and drain water that is in the washer. The alarm is canceled by opening the door. See *Chapter 7: Troubleshooting – Washer leaks.*

### LIQ DETER

This displays only if the optional liquid detergent dispenser is installed and the level of the detergent in the dispense container is low. The alarm is disabled if the detergent dispense volume is set to “0.” The alarm is canceled when detergent is added to the container.

### DRAIN FAIL

This displays if water fails to completely drain during a normal drain operation. Operation of the washer is canceled. The alarm is canceled by opening the door.

### LOW TEMP

This displays if the water doesn't reach the set point temperature within 135 minutes. Operation of the washer continues. The alarm cancels when the door is opened.

---

# CHAPTER 5

## MAINTAINING YOUR WASHER

Under normal operation, your Glassware Washer will require little routine maintenance to keep it functioning properly.

### **Maintaining the Stainless Steel Interior**

The interior of your Glassware Washer is type 304 stainless steel. Never use metal scouring pads on the stainless steel interior. Metal scouring pads can scratch or leave metal chips that will rust.

To maintain the appearance and quality of the stainless steel interior, perform the following tasks as needed:

- Wipe the outer edges on the inside door panel to remove particles and residue that occur during loading.
- Use a sponge sprinkled with non-abrasive stainless steel cleaner to gently wipe the washer tank and door to remove surface discoloration caused by exposure to halogenated compounds or steel parts. In some cases, the finish may be slightly dulled by this action.

If your water has a high mineral content, film or spots may develop on the washer interior. If the procedures listed above do not eliminate the film or spots, you may run your washer through the special cycle below.

### **To remove film or spots from the washer interior:**

1. Remove all glassware from the washer.
2. Choose the Glass Program, Steam OFF, Pure Rinse OFF, and set the drying time to zero.

3. After the first fill is complete, open the washer and carefully pour one cup of bleach into the bottom of the tank.
4. Close the washer door and allow the unit to wash and drain.
5. When the washer fills the second time, open the washer and add two cups of 4-8% acetic acid (or white vinegar) to the bottom of the tank.
6. Close the washer door and allow the unit to wash and drain without further interruption.


### Maintaining the Air Gap in Drain

If the installation included an air gap (recommended) it requires periodic cleaning. The air gap is not part of the washer; therefore, it is not covered by the washer warranty. The air gap protects the washer against water backing up in the event of a clogged drain. Check the air gap for build-up any time your washer is not draining well.

Most types of air gaps are easy to clean. Simply follow the manufacturer's directions for maintaining your air gap.

### Maintaining the Water Fill Valve

The filter screen of the water fill valve, illustrated below, may require periodic cleaning to remove trapped particles. A low water fill occurs if the blockage is not removed.

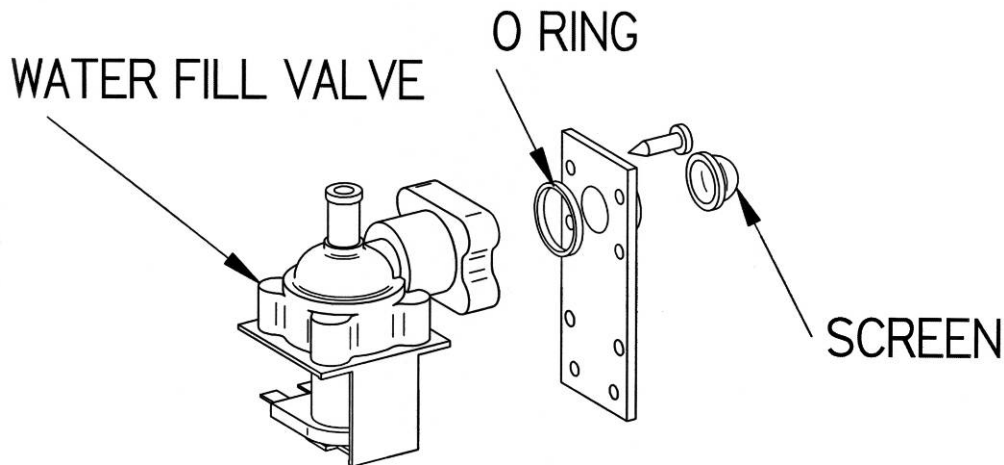

#### **To clean the water fill valve:**

1. Turn off the water and electrical power to the washer. Unplug the Mobile model; turn off the circuit breaker that services the Undercounter or Freestanding model.
2. Remove the rear panel of FreeStanding models.
3. Disconnect the water supply tube that is connected to the valve.
4. Remove the valve from the base.


5. Remove the four screws from the valve plate and separate the valve from the inlet connection.
6. Remove the valve filter screen and clean it under running water.
7. Replace the screen and reassemble the valve connection. Be certain the O-ring gasket is positioned in the groove in the fill valve body.

## Replacing the Washer Light Bulb

If your Glassware Washer is equipped with a window, you will need to periodically replace the light bulb. Refer to the figure below and follow the instructions below to replace the washer light bulb.

### To replace the washer light bulb:

1. Turn off electrical power to the Glassware Washer.
2. Remove the lower and middle panels as described in *Chapter 3: Getting Started, Connecting the Power Supply*.
3. Disconnect the springs on each side that connect the pivot brackets to the washer base. Disconnect the hook on the washer base side.
4. Remove the two screws on each side of the outer door panel. Lower the panel and remove it.
5. Remove the bulb by pressing in on the bulb and rotating the bulb counterclockwise.
6. Insert the new bulb and rotate clockwise into place.
7. Replace the outer door panel.

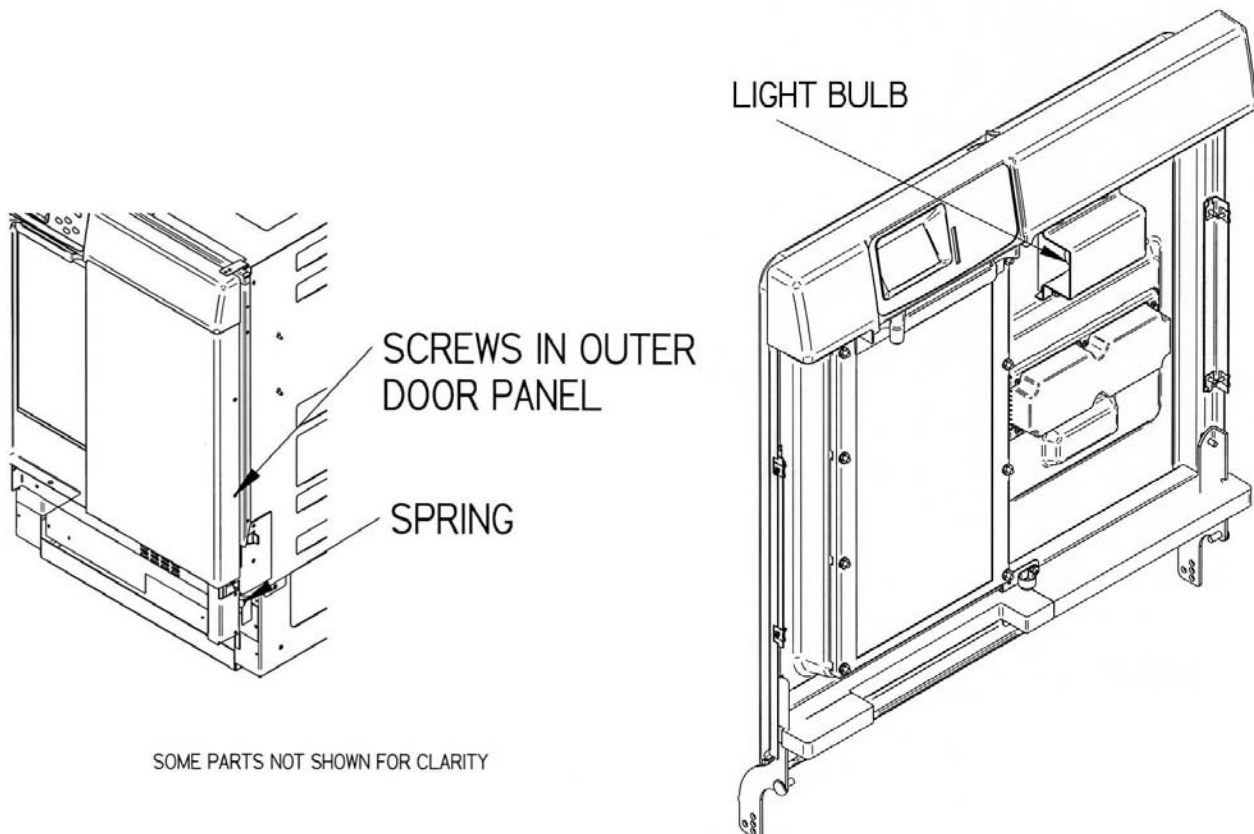


## Removing the Washer Top, Back, and Side Panels

There may be times when you need to remove the washer top, back, and side panels to clean, maintain, or repair your washer. Refer to the figure below and follow the instructions to remove the top, back, and side panels of your washer.

### **To remove the washer top, back, and side panels:**

1. Turn off electrical power to the Glassware Washer.
2. Remove all of the screws from the back panel of the washer and remove the back panel.
3. Remove the 2 screws at the kick panel.
4. Remove 2 screws at the rear that hold the sides to the top.
5. Remove the two screws from each side of the washer at the front edge.
6. Remove the side panels.
7. Remove the 2 screws that hold the top.
8. Remove the top.

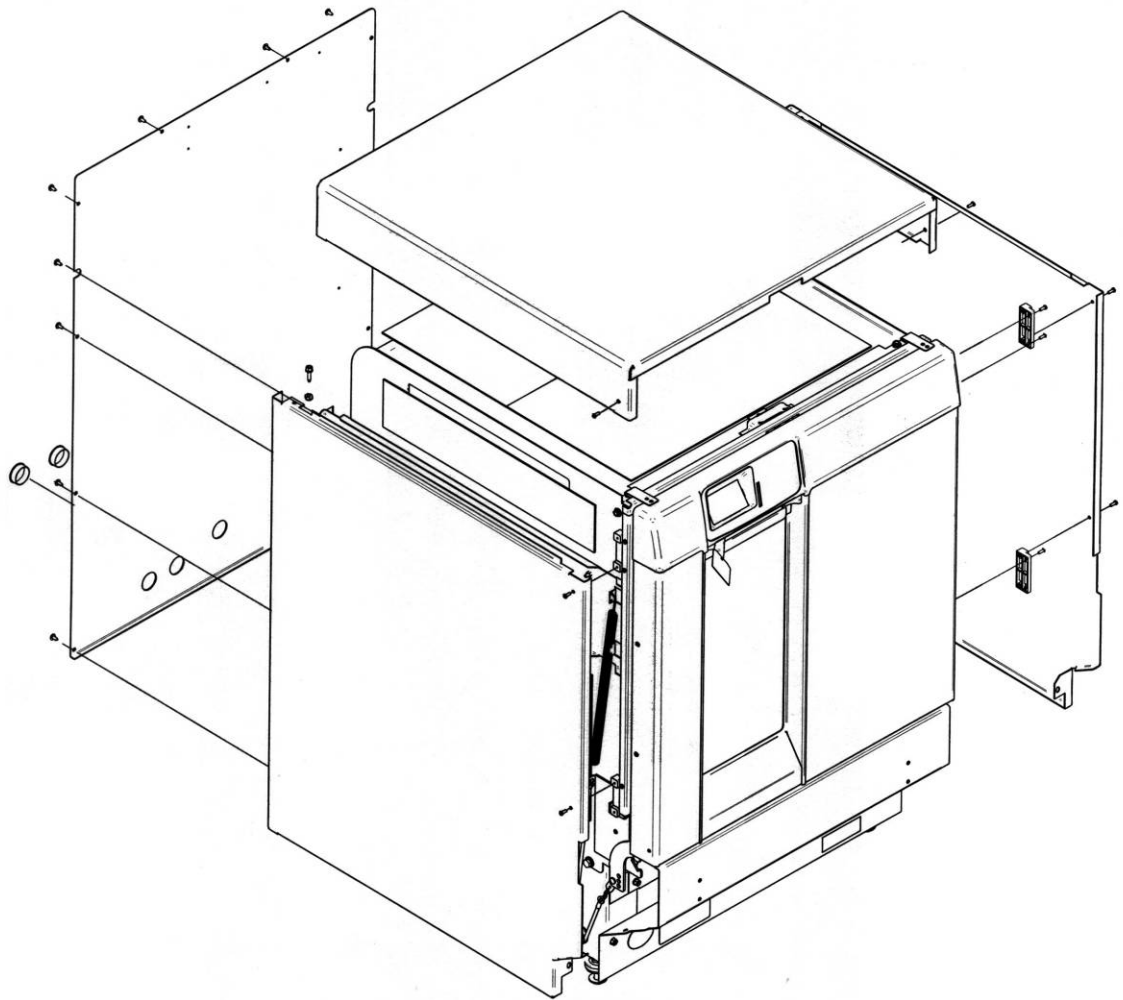


## Cleaning the Sump Filter Screen

If debris accumulates on the filter screen that is located in the washer sump, you can remove and clean the screen. To clean the sump filter screen, refer to the figure below and follow the instructions below.

### To clean the sump filter screen:

1. Open the washer door and remove the bottom rack.
2. Remove the screw that retains the filter screen.
3. Carefully lift the screen and pull forward, using care not to drop debris into the sump.

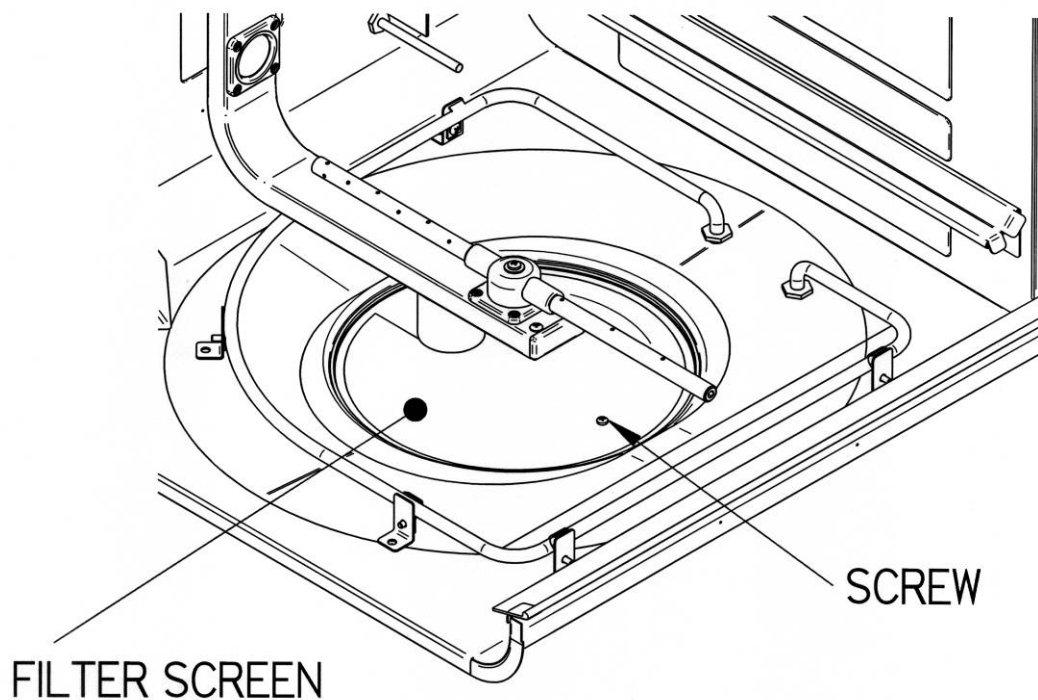

---


# CHAPTER 6

## MODIFYING YOUR WASHER

The configuration of your washer may be changed to accommodate your needs. For example, you may want to install pipet inserts, install an optional top rack, or install an optional liquid detergent dispense system or Cool Drain Water Kit. Refer to *Appendix A: Glassware Washer Accessories* for the accessories that are available for your Glassware Washer.

### Adjusting the Height of the Upper Rack

The height of the upper rack may be varied to suit the heights of glassware that will be washed in the upper and lower racks.

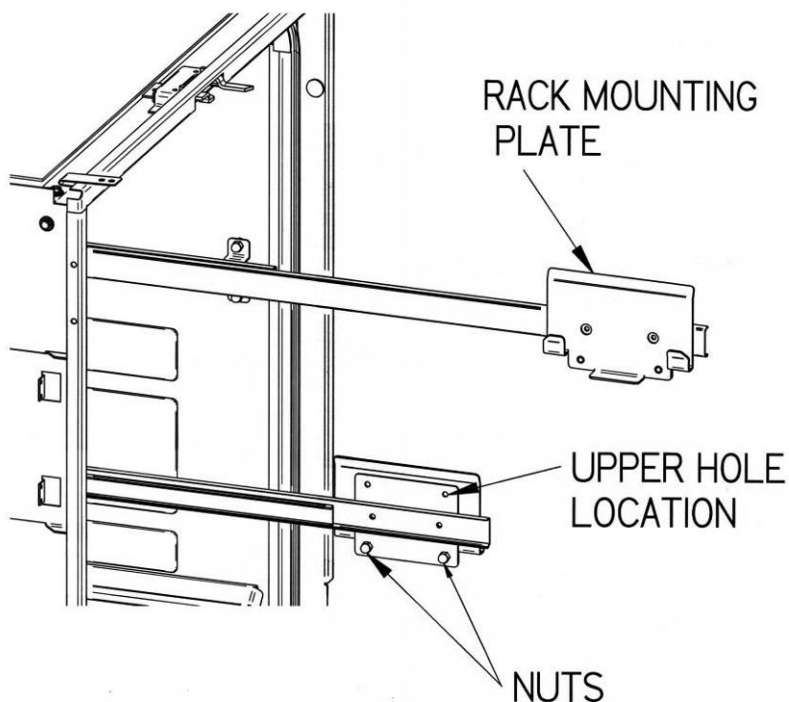



**To adjust the height of the upper rack:**

1. Remove the upper rack as described in *Chapter 4: Using Your Washer*.
2. Remove the 2 nuts that retain the rack mounting plate to the slide.
3. Reposition the rack mounting plate to the slide using the other set of holes.  
Replace the 2 nuts.
4. Repeat the procedure on the opposite slide.

**Spindle Rack Options**


If you purchased a FlaskScrubber Glassware Washer, you can modify the configuration of your washer to better suit your needs. Optional accessories include two Pipet Inserts, and an Upper Spindle Rack. The Pipet Inserts are easily installed in place of one or more spindles in the washer and are designed to accommodate 1 ml to 50 ml pipets.

**Installing Optional Pipet Inserts in a FlaskScrubber**

Two sizes of Pipet Inserts are available. The 8-place Pipet Insert, shown below accommodates 1 ml to 10 ml pipets. Multiple Pipet Inserts may be installed, if desired. Refer to the figure and follow the instructions below to install an 8-place Pipet Insert.

**To install an 8-place Pipet Insert:**


1. Insert the tool supplied with the Spindle Rack or a screwdriver into the top of a spindle and turn counterclockwise to remove the spindle.
2. In the position previously occupied by the spindle, thread into place the short spindle you received with the Pipet Insert.
3. Push the Pipet Insert over the short spindle.
4. If spindles other than the one used for the short spindle are removed, the holes must be plugged with the plastic plugs supplied with the spindle rack.

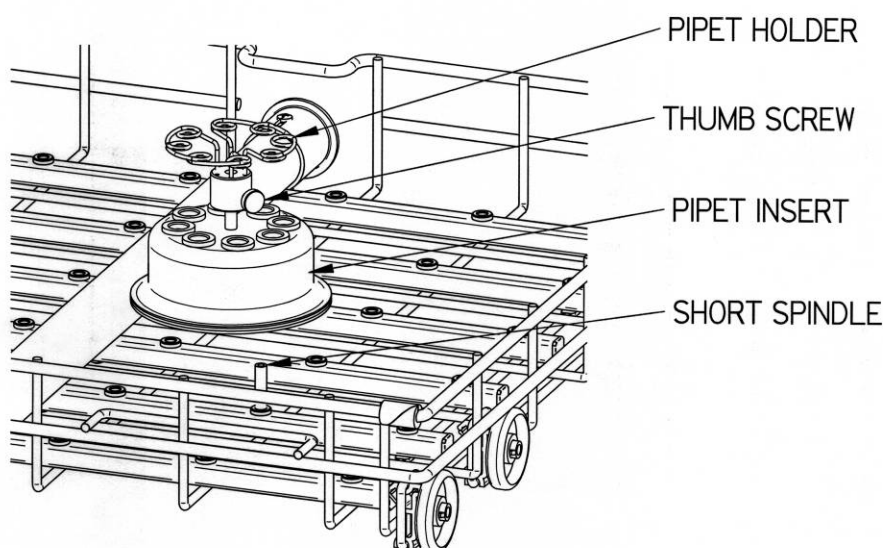

The 16-place Pipet Insert, shown below accommodates 1 ml to 50 ml pipets. Two of these inserts may be installed, if desired. Refer to the figure and follow the instructions below to install a 16-place Pipet Insert.

### **To install a 16-place Pipet Insert:**

1. Use the tool supplied with the Spindle Rack or a screwdriver to remove the six spindles on either the front or rear of the spindle rack. Remove others if necessary.
2. Replace the left and right spindles with the two short spindles that you received with the Pipet Insert.
3. Replace all other spindles with the plastic plugs that you received with your washer.
4. With the support facing the front of the spindle rack, slide the Pipet Insert over the spindles until it completely seats.
5. Push pipets into the 16 receptacles and rest them on the support.

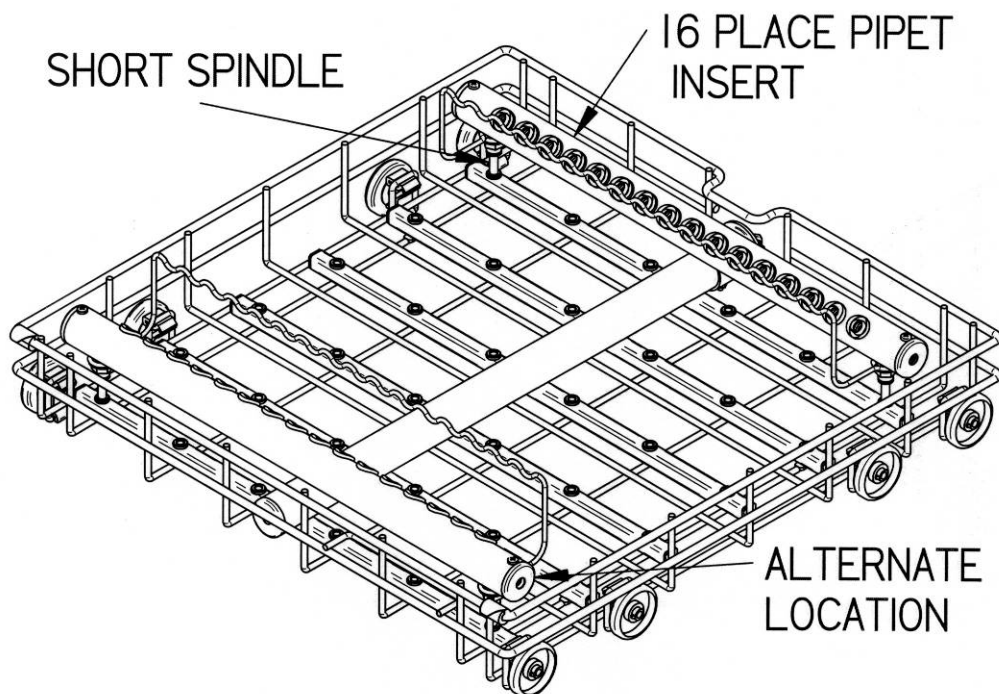

**Note:** Lower spindle racks are shipped with 34 long spindles and 2 shorter spindles. If spindles are removed and then replaced, the shorter spindles should be located on the front row in the second and third positions from the right side. This will allow proper function of the detergent dispenser.


## Installing the Optional Upper Spindle Rack

The Upper Spindle Rack expands the capability of the FlaskScrubber by providing 30 additional spindles for washing narrow-neck glassware. The rack may be customized by removing spindles or by installing Multi-Pin Inserts which hold beakers or other wide mouth glassware.


### To install the optional Upper Spindle Rack:

Install the upper rails and the upper spindle rack. See the figure below.

1. Remove the acorn nuts and lockwashers from both upper sides using a 3/8" wrench. There are 4 on each side.

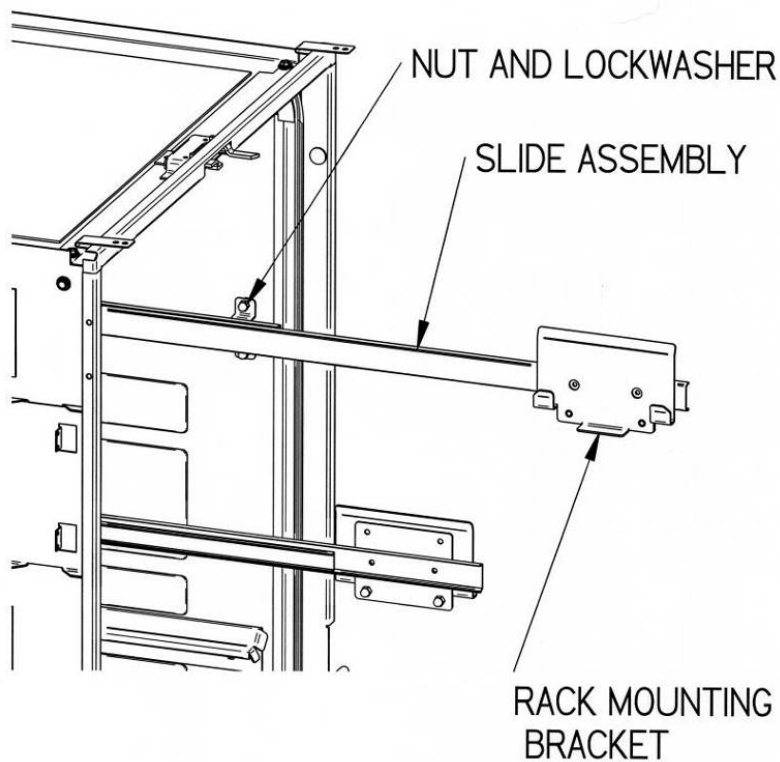

2. Position brackets on slides over the studs on the sides. The position of the rack mounting plates must be as shown. Secure the slide assemblies using lockwashers and acorn nuts removed in Step 1. It may be necessary to synchronize the travel of the two slides. Pull each slide out firmly to their maximum travel.

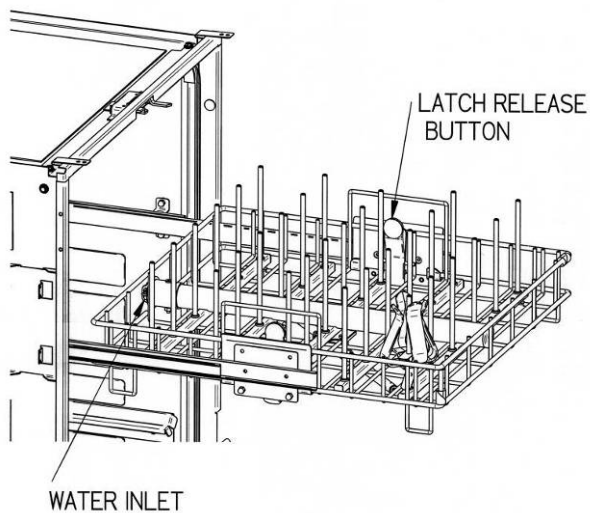

3. Install the Upper Spindle Rack by lowering the rack onto the rack mounting plates. The water inlet must have its inlet toward the rear. Make sure the latch on each side secures the rack to the rack mounting plate.


When the Upper Spindle Rack is in place, it limits the height of the glassware that can be washed. If very tall glassware or long pipets are to be washed, remove the Upper Spindle Rack.

### Installing Optional Inserts into the Spindle Rack

A Multi-Pin insert is available that can be placed directly on the FlaskScrubber upper or lower spindle rack. It can hold beakers or wide mouth glassware. The insert can be positioned either on the left side or right side.

#### **To install an insert:**

1. Remove all glassware, spring retainers and glassware holders.
2. Lower the insert down over the spindles. If necessary, spindles may be removed using the tool supplied with the Spindle Rack or a screwdriver in the slot on the top of the spindles. Be sure to replace any removed spindle with a plug supplied.

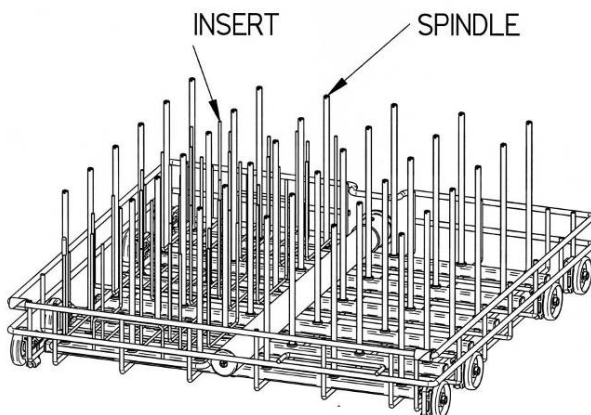

## Installing Small Spindles

Ten small spindles are included with each Spindle Rack. Since the large volume of water which flows through the standard spindles may be great enough to force the glassware off the spindle, small spindles should be used when small pieces of glassware will be washed.

### **To install small spindles:**


1. Remove the standard spindle by inserting the tool supplied with the Spindle Rack or a screwdriver into the slot on the top and turning counterclockwise.
2. Replace the standard spindle with a small spindle.

## Installing the Optional Cool Drain Water Kit

An optional Cool Drain Water Kit is available to reduce the water temperature of the drain water in installations where hot drain water is not permitted. It may be installed on Undercounter or Freestanding models. A supply of cold water must be available. It is recommended that the cold water temperature is colder than 54°F (12°C). Warmer water temperature will cause the drain temperature to be higher. The cold water supply must be capable of a flow rate of 3.5 GPM (13.2 LPM) and a pressure of 20-120 PSI (138-827 kPa). Hot drain water will be mixed with cold water during drain operations. To install the optional Cool Drain Water Kit, refer to the figure below and the following instructions.

### **To install the optional Cool Drain Water Kit:**

1. Disconnect power to the washer.
2. Remove the back panel of the washer if it is a Freestanding model or the cover panel of an Undercounter model.
3. Attach the water valve assembly to the rear flange of the base using the screws provided.
4. Connect the wires coming from the water valve assembly to the mating connector in the wiring harness of the washer.
5. Connect the cold water supply to the water valve assembly. Do not allow the tube to kink when the washer is placed into its operational location. If necessary, replace the straight fitting on the cool water valve with the supplied elbow.
6. Replace the back panel if it is a Freestanding model or the cover panel if it is an Undercounter model.
7. Attach the hose from the valve to a Twin Inlet High Flow Air-gap suitable for dishwasher applications. This may be obtained from a plumbing supply store.
8. Attach the drain hose from the Glassware Washer to the other inlet of the Air-gap.
9. Attach the Air-gap to the building drain.
10. Turn on the electrical power.

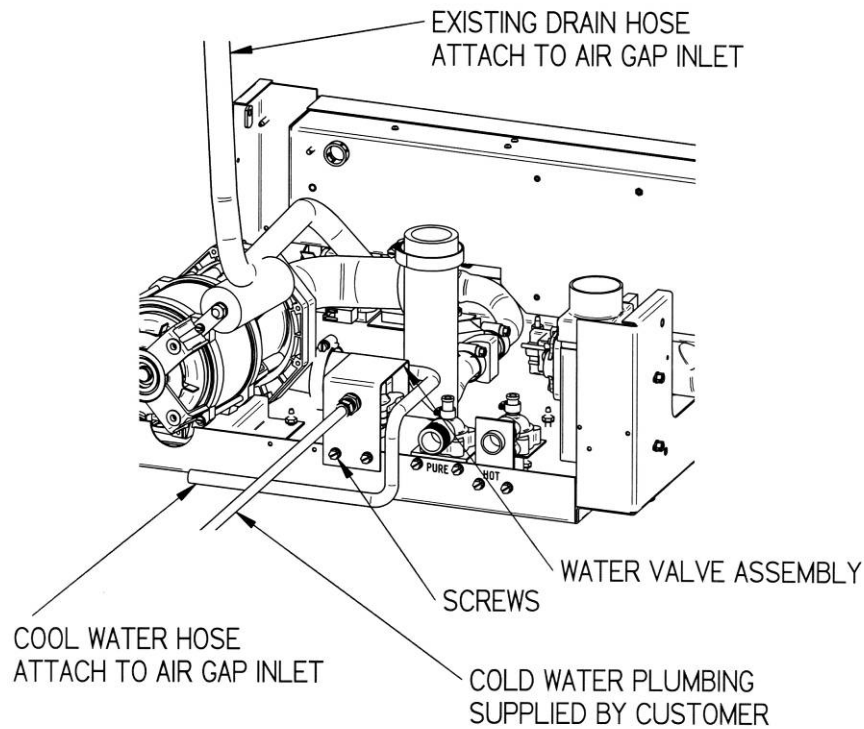

TANK AND OTHER PARTS NOT SHOWN FOR CLARITY


## Installing the Optional Liquid Detergent Dispenser


An optional liquid detergent dispenser is available for installation on SteamScrubbers and FlaskScrubbers. This permits dispensing liquid detergent from a bulk 1 gallon (3.8 L) container eliminating the need to fill the detergent cup with detergent each time the washer is used. To install the liquid detergent dispenser, refer to the figures below and the following instructions.

### **To install the liquid detergent dispenser components:**

1. Disconnect power to the washer.
2. Remove the back panel if the washer is a Freestanding model. Loosen the insulation blanket to provide access to the back of the washer.
3. Attach the dispense pump to the rear flange of the base using the screws and nuts provided.
4. Attach the tube from the detergent container to the inlet of the pump.
5. Loosen the hose clamp that is around the large hose on the bottom of the vacuum break. Do not remove the hose. Slip the stainless steel tube that is on the end of the clear tube coming from the detergent dispense pump into the hose on the vacuum break. Tighten the hose clamp. Secure the provided tie wrap around all 3 tubes on the top of the vacuum break being careful not to collapse any tube.
6. Locate the detergent container at a convenient location where it will not be a trip hazard.

7. Remove the lower panel and the electrical junction box cover as described in *Chapter 3: Getting Started*.
8. Plug the supplied harness into terminal J6 on the Relay PCB and route the other end to the left rear corner of the washer base.
9. Plug the wire from the detergent container to the mating connector in the new harness in the base of the washer.
10. Plug the connectors on the pump to the mating connectors in the new harness in the base of the washer.
11. Reinstall the junction box cover and lower panel.
12. Reinstall the back panel if the washer is a Freestanding model.
13. Turn on the electrical power.

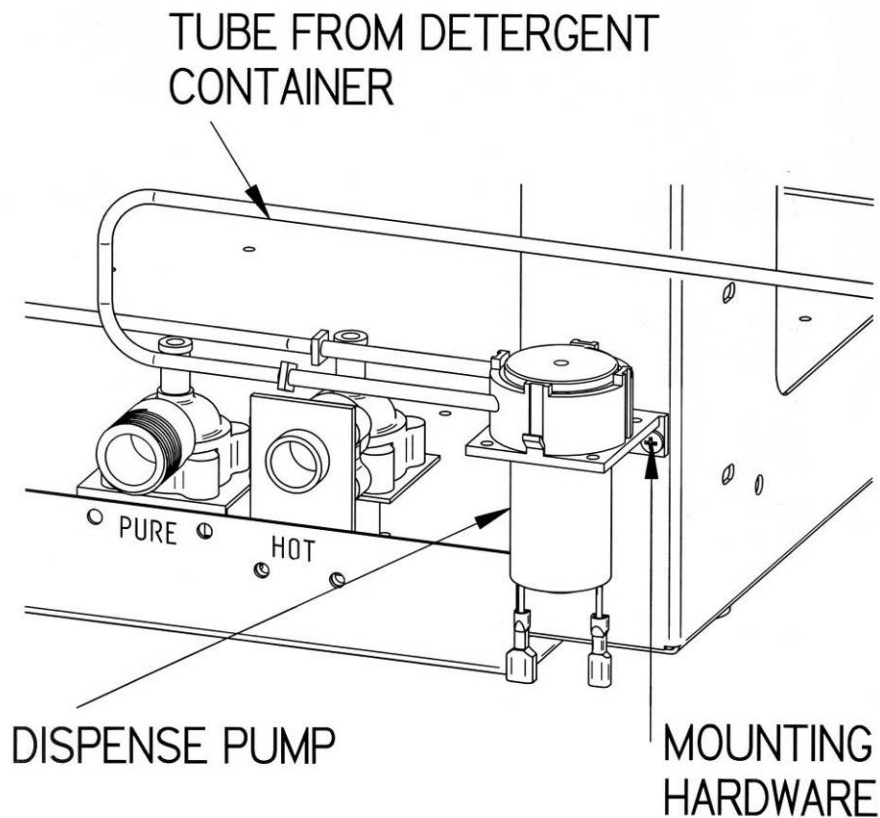

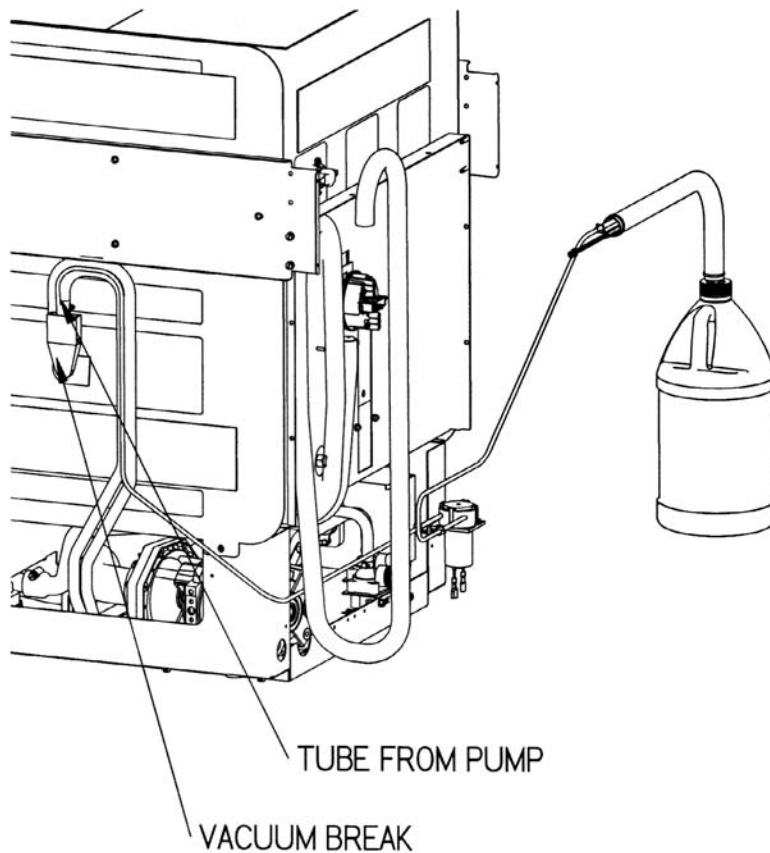

**Enable the liquid detergent dispenser operation:**

1. Enter the washer set up screen on the display by holding the ▼ button on the control panel and closing the door and sliding the door latch to the right.
2. Press the ▼ button until the arrow on the left points to LIQ DETERGENT OPT.
3. Press the ► or ◀ button to activate this accessory.
4. Press RUN to save all settings.

**Using the liquid detergent dispenser:**

1. Remove the cap from the container and remove the tube and level switch assembly.
2. Fill the container with detergent or use a new bottle of Labconco detergent and replace the tube and switch assembly and the cap.
3. There are two methods which can be followed for filling the system with detergent.
  - a. Run a wash cycle with no glassware in the washer. Set the detergent dispense rate to 30 ml.
  - b. Run a diagnostics test. When the display window shows DETERGENT PUMP, allow the pump to run for 30 seconds.
4. Program the quantity of detergent that will be dispensed by entering the programming screen as described in *Chapter 4: Using Your Washer*.



---

# CHAPTER 7

## TROUBLESHOOTING

Refer to the following table if your Glassware Washer fails to operate properly. If the suggested corrective actions do not solve your problem, contact Labconco Product Service for additional assistance.

| PROBLEM                                     | CAUSE   | CORRECTIVE ACTION   |
|---|---|---|
| <b>Detergent cup fails to open</b>          | Malfunction of cup                                | Replace cup.  |
|   | Malfunction of relay                              | Check electrical continuity or replace PCB.   |
|   | Glassware interfering with cup door               | Move glassware.   |
| <b>Noises during operation</b>              | Glassware rattling                                | Load glassware properly. Do not allow glassware to touch.   |
|   | Door panel vibrating                              | Tighten panel screws.   |
|   | Wash arms hitting                                 | Check glassware loading.  |
|   | Water inlet line makes occasional hammering sound | Check with plumber.   |
|   | Purified water pump is activated                  | No corrective action needed. This is normal for about 5 minutes while pure water enters the washer. |
| <b>Door drops too far (below door stop)</b> | Bent door hinge                                   | Replace door hinge.   |

| PROBLEM                      | CAUSE   | CORRECTIVE ACTION  |
|------------------------------|---|--|
| <b>Poor drying</b>           | Dry temperature too low   | Program warmer dry temperature.  |
|                              | Flask, beaker, or labware with concave bottoms not loaded correctly | Tilt these articles when loading so that water drains off.   |
|                              | Recondensation  | If clean labware is left in the washer, recondensation may occur. Open latch after completion of cycle to allow moisture-laden air to escape.  |
|                              | Heater(s) disconnected or malfunctioned                             | Reconnect or replace heater.   |
|                              | Blower disconnected or malfunctioned                                | Reconnect or replace blower.   |
|                              | Dry time too short  | Program longer time.   |
|                              | Temperature sensor out of calibration                               | Verify displayed temperature by attaching a thermocouple to the temperature sensor. Route leads outside of the washer. Close and latch the door while holding the ◀ button. Call Labconco Product Services if the displayed temperature differs from the test temperature by more than $\pm 4^{\circ}\text{C}$ . |
| <b>Poor washing results</b>  | Water is not hot enough   | Check Wash 2 water temperature. Must be at least 120°F (49°C). Recommended temperature is 150°F (66°C).  |
|                              | Improper loading  | Check washer loading instructions in <i>Chapter 4</i> in this manual.  |
|                              | Detergent too old or amount wrong                                   | Replace detergent or change quantity used.   |
|                              | Detergent caked in cup  | Clean cup. Replace with fresh detergent. Check water inlet temperature. Must be at least 120°F (49°C).   |
|                              | Insufficient wash time  | Increase wash cycle time.  |
|                              | Insufficient water  | Check to be sure the water shut-off valve is fully open. Check to be sure the water pressure is between 20 and 120 psi (138-827 kPa) at a flow rate of 1.25 (4.7 Liters/min) gallons per minute. Check water level switch.   |
| <b>Not draining properly</b> | Clogged filter  | Clean sump filter.   |
|                              | Drain line clogged or kinked  | Disconnect drain line and clean or remove kink.  |

| <b>PROBLEM</b>                               | <b>CAUSE</b>                                  | <b>CORRECTIVE ACTION</b>   |
|--|---|--|
| <b>Purified water does not fill</b>          | Crimp in supply hose                          | Refer to installation instructions for proper purified water piping installation.                                      |
|  | Motor wires disconnected                      | Reconnect wires.   |
| <b>Tank and inner door have surface rust</b> | Collects normally                             | Clean with stainless steel cleaner and soft cloth.   |
|  | Glassware is contaminated with mineral acids. | Pre-rinse glassware. Clean with stainless steel cleaner and soft cloth.  |
| <b>Washer does not run</b>                   | No power to washer                            | Check fuse or circuit breaker, or door switch.   |
|  | Door latch is open                            | Latch washer door by pushing door in.  |
| <b>Washer door opens and closes badly</b>    | Cabinet opening is not square                 | Correct opening with leveling feet.  |
|  | Door hitting counter-top mounting screws      | Adjust mounting screws.  |
| <b>Washer leaks</b>                          | Washer is not level                           | Adjust leveling feet.  |
|  | Door not sealing                              | Adjust door latch.   |
|  | Tank gasket loose                             | Check gasket and correct.  |
|  | Loose connection at hose clamps               | Tighten hose clamp.  |
|  | Loose connection at fill valve                | Check fill valve and correct.  |
|  | Improper detergent                            | Use non-foaming detergent. Use only proper amount of detergent.  |
|  | Obstruction in drain line                     | Remove obstruction.  |
|  | Washer overfilling                            | Check operation of fill switch.  |
| <b>Washer leaks around door seal</b>         | Improper installation                         | Check cabinet opening to see if it is aligned properly with the door. Washer must be level. Adjust feet to level unit. |
|  | Unit not level                                | Adjust feet to level unit.   |
|  | Door opened during operation                  | Allow 4-5 seconds before opening door after unlatching.  |

| PROBLEM  | CAUSE  | CORRECTIVE ACTION  |
|--|--|--|
| <b>Washer will not drain</b>   | Clogged drain  | Clean out drain.   |
|  | Blockage of air gap  | Clean the air gap by lifting the cover, unscrewing the cap, and removing any accumulated material. See <i>Chapter 5</i> .  |
|  | Kink in drain hose   | Check for a kink in the drain hose.  |
|  | Filter blocked   | Check filter in washer and clean as needed.  |
|  | Drain valve or pump not operating  | Check operation.   |
| <b>Washer will not fill</b>  | Fill valve clogged   | Clean fill valve.  |
|  | Water supply turned off  | Turn on water.   |
|  | Defective fill switch  | Check operation.   |
| <b>Washer overfills</b>  | Defective fill switch  | Check operation.   |
| <b>Selection buttons on control panel inoperable or controls malfunction</b> | Microprocessor memory error  | Perform "Factory Reset." Open latch on door. Press ▼ button and simultaneously close door latch. Release ▼ button and display will show the set up screen. Press ▼ until FACTORY RESET is highlighted. Press ► or ◀ until "Yes" is displayed. Press RUN. |
|  | Key pad defective  | Replace key pad.   |
| <b>Drain Fail Alarm</b>  | Drain Valve inoperable   | Listen for click when washer should drain.   |
|  |  | Check for mechanical restriction of actuator arm on top of valve.  |
|  |  | Check for electrical power to valve.   |
|  |  | Check for damage to valve coil.  |
|  | Drain Pump inoperable  | Check if pump fan turns when washer should drain.  |
|  |  | Check for mechanical restriction of fan.   |
|  |  | Check for electrical power to pump.  |
| Drain hose not routed properly   | Hose from drain pump must slope upward and have no low spots or horizontal runs under washer tank. |  |
|  | Make sure hose is attached to the back of the washer and has a high loop.                          |  |

| PROBLEM                 | CAUSE                                  | CORRECTIVE ACTION  |
|-------------------------|--|--|
|                         | Drain hose not routed properly (cont.) | Hose must be attached to building plumbing pipe $\frac{3}{4}$ inch ID or larger.<br><br>Hose from washer must not be attached to another hose.<br><br>Check for kinks in hose.<br><br>Drain hose must not be elevated more than 30 inches. |
|                         | Pump lost its prime                    | Remove ALL water from inside drain hose. Pump will prime when water drains from washer.  |
|                         | Building plumbing defective            | Check for blocked air gap.<br><br>Check for clogged drain.   |
|                         | Filter screen clogged                  | Clean screen in bottom of washer.  |
|                         | Defective fill level switch            | Contact Labconco Product Service.  |
|                         | Defective controller                   | Check output from PCB to drain valve and pump.<br><br>Replace PCB. Contact Labconco Product Service.   |
| <b>Water High Alarm</b> | Holes in filter screen too small       | Contact Labconco Product Service.  |
|                         | Defective hose to fill switch          | Check for splits.<br><br>Check for loose clamps.   |
|                         | Defective fill level switch            | Recalibrate switch. Contact Labconco Product Service.<br><br>Replace switch.   |
|                         | Holes in filter screen too small       | Contact Labconco Product Service.  |
| <b>Water Low Alarm</b>  | Improper water supply                  | Make sure water is turned on.<br><br>Tap water pressure must be at least 20 psi (138 kPa) and flow must be at least 1.25 gallons (4.7 liters)/minute.<br><br>DI water supply must have flow of at least 0.9 gallons (3.4 liter)/minute.    |



| PROBLEM               | CAUSE  | CORRECTIVE ACTION   |
|-----------------------|--|---|
|                       | Improper water supply (cont.)                              | At least 3.4 gallons (13 liters) must be available for each DI rinse.   |
|                       | Defective fill valves                                      | Clean screens inside the inlet fittings of the water valves.<br><br>Check voltage to valves.<br><br>Replace valve if it does not operate. |
|                       | Restricted flow  | Check for kinks in water hoses in base of washer to water inlet fitting on side of tank.  |
|                       | Defective DI water pump                                    | Check voltage at pump.<br><br>Listen for pump operation. Replace if necessary.  |
|                       | Defective hose to fill switch                              | Check for splits.<br><br>Check for loose clamps.  |
|                       | Defective fill level switch                                | Recalibrate switch. Contact Labconco Product Service.<br><br>Replace switch.  |
|                       | Defective controller                                       | Check output from PCB to both valves and DI pump.<br><br>Replace PCB. Contact Labconco Product Service.                                   |
|                       | Holes in filter screen too small                           | Contact Labconco Product Service.   |
|                       | Defective grommet where hose connects to fill level switch | Replace grommet.  |
| <b>Low Temp Alarm</b> | Heater inoperative   | Check electrical connections.<br><br>Check voltage to heater.<br><br>Check resistance of heater. If open circuit, replace heater.         |
|                       | Defective controller                                       | Check output from PCB to heater.<br><br>Replace PCB. Contact Labconco Product Service.  |

| <b>PROBLEM</b> | <b>CAUSE</b>                              | <b>CORRECTIVE ACTION</b>  |
|----------------|---|---|
|                | Defective temperature probe               | Check electrical connections.<br><br>Contact Labconco Product Service. Replace if necessary.    |
|                | Defective high temperature cut out switch | Check resistance of safety cut out switch. If open circuit at room temperature, replace switch. |




# APPENDIX A

## GLASSWARE WASHER ACCESSORIES




### FlaskScrubber Accessories




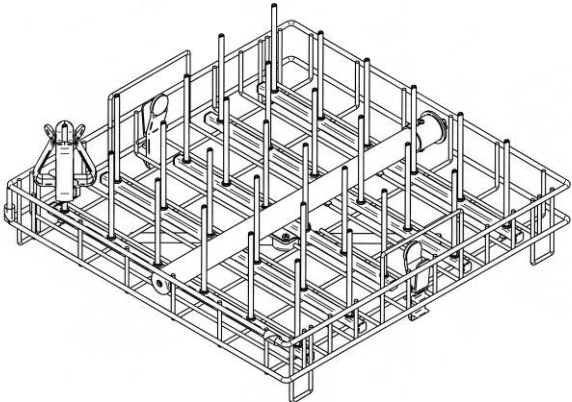
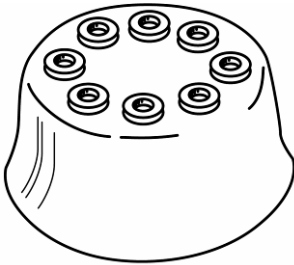
| PART #  | DESCRIPTION   |   |
|---------|---|---|
| 4595600 | <b>Upper Spindle Rack – Stainless Steel.</b> Contains one upper spindle rack with 30 spindles, 30 glassware holders and 2 slides. |    |
| 4595800 | Same as above except without slides.  |   |
| 4494200 | <b>8-Place Pipet Washer – Stainless Steel.</b> Holds 1 ml to 10 ml pipets.  |  |

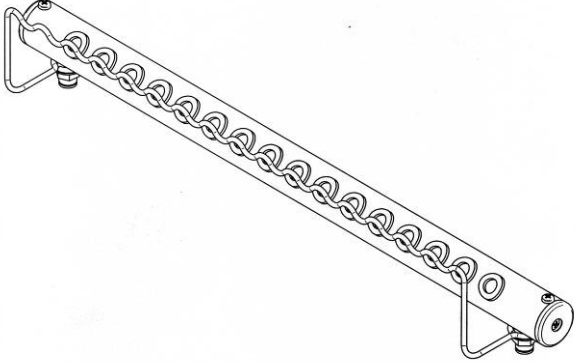
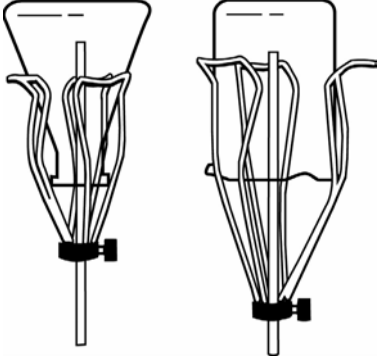
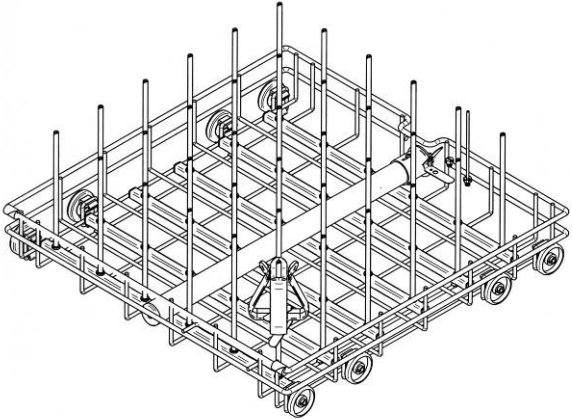





| PART #                        | DESCRIPTION   |  |
|-------------------------------|---|--|
| 4595200                       | <p><b>16-Place Pipet Washer – Stainless Steel.</b> Holds 1 ml to 50 ml pipets.</p>  |    |
| <p>4424600</p> <p>4424800</p> | <p><b>Glassware Holder – Small.</b> Grips wide-mouth glassware over spindles. Vinyl coated stainless steel wires.</p> <p><b>Glassware Holder – Large.</b> Grips wide-mouth glassware over spindles. Vinyl coated stainless steel wires.</p> |    |
| 4596000                       | <p><b>Small Spindles with Clips.</b> Replaces larger spindles when washing small glassware.</p>   |  |
| 4595700                       | <p><b>Lower Spindle Rack – Stainless Steel.</b> Contains 36 spindles and 36 glassware holders. Included with FlaskScrubbers.</p>  |  |



## FlaskScrubber or SteamScrubber Accessories
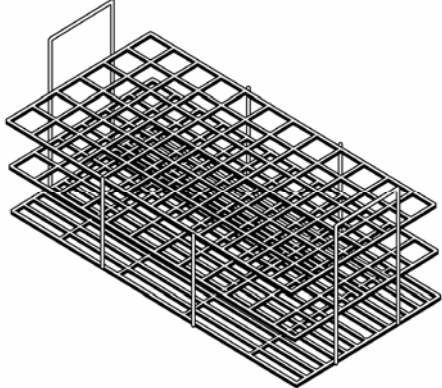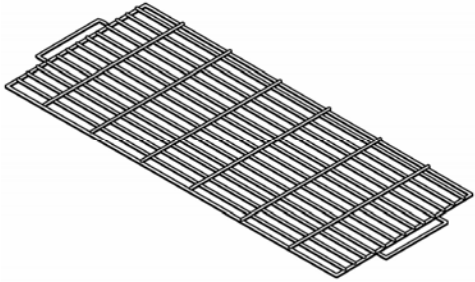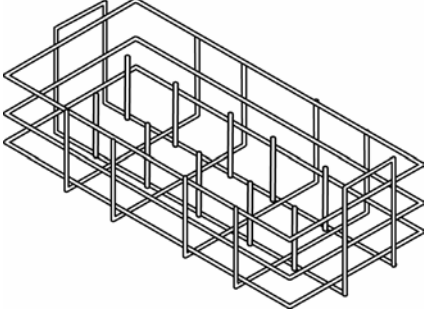
| PART #  | DESCRIPTION  |  |
|---------|--|--|
| 4402001 | <p><b>Culture Tube/Test Tube Insert – Stainless Steel.</b><br/>Insert for 15 mm to 18 mm tubes.</p>  |    |
| 4402101 | <p><b>Culture Tube/Test Tube Insert – Stainless Steel.</b><br/>Insert for 20 mm to 25 mm tubes.</p>  |  |
| 4401301 | <p><b>Culture tube/Test Tube Insert – Stainless Steel.</b><br/>Insert for 10 mm to 12 mm tubes. Includes retainer top 4401401.</p>                                   |  |
| 4401401 | <p><b>Retainer Top for Culture Tube Insert – Stainless Steel.</b> For use with all culture tube inserts to keep lightweight tubes in place under water pressure.</p> |   |
| 4401501 | <p><b>10-Pin Insert – Stainless Steel.</b> For use with beakers, Erlenmeyer flasks, and other wide-mouth glassware.</p>  |  |
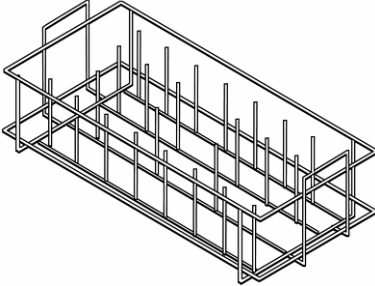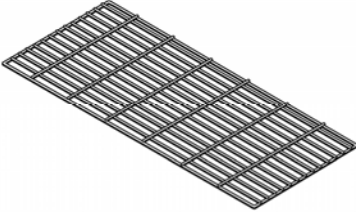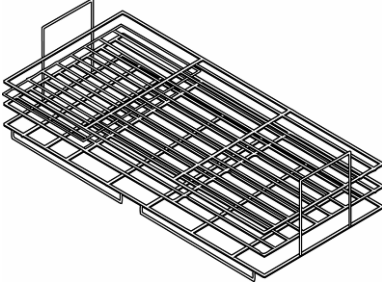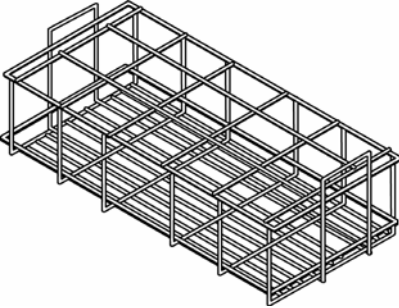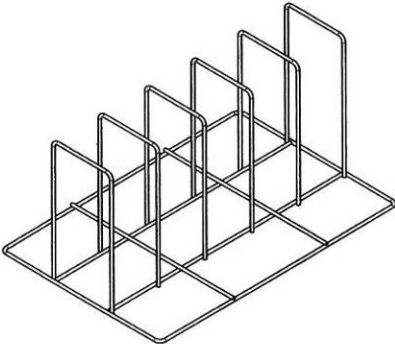
| PART #  | DESCRIPTION   |  |
|---------|---|--|
| 4401801 | <p><b>32-Pin Insert – Stainless Steel.</b> For use with beakers, Erlenmeyer flasks, and other wide-mouth glassware.</p>   |    |
| 4401901 | <p><b>Retainer Top – Stainless Steel.</b> To keep lightweight flasks in place under water pressure. For use with 4401801 or 4401501.</p>                              |    |
| 4589701 | <p><b>Petri Dish Insert – Stainless Steel.</b> For plates, petri dishes, and watch glasses.</p>   |   |
| 4589201 | <p><b>BOD Rack/Beaker/ Erlenmeyer Flask Insert – Stainless Steel.</b> Rack secures glassware used in oxygen-demand determination, beakers, and Erlenmeyer flasks.</p> |  |
| 4598401 | <p><b>Tray Holder.</b> For trays up to 19 x 19 x 2.5 deep (48 x 48 x 6 cm deep)</p>   |  |
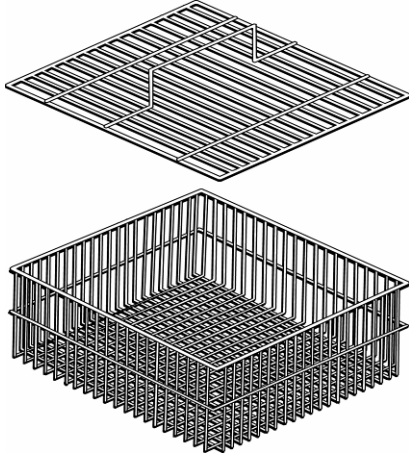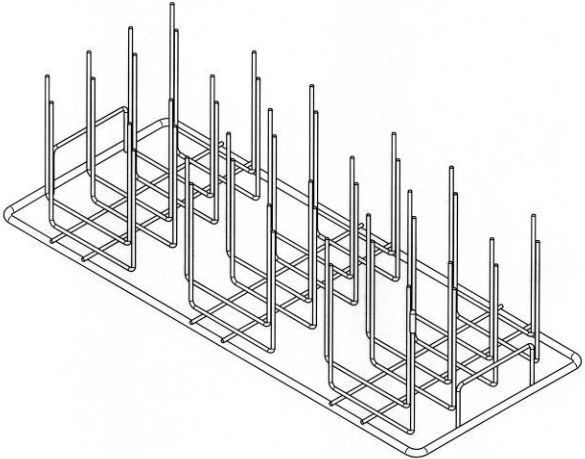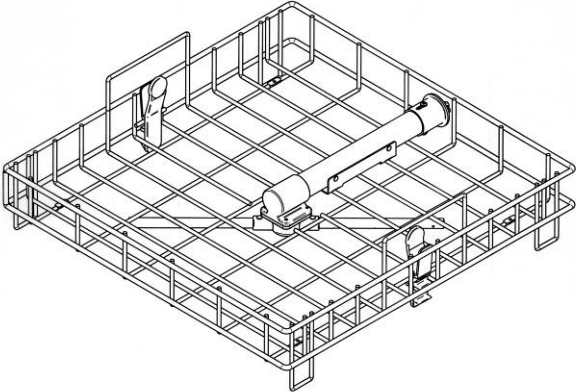
| PART #                 | DESCRIPTION   |   |
|------------------------|---|---|
| 4402201                | <p><b>Utensil Basket – Stainless Steel.</b> For miscellaneous labware such as stoppers, spatulas, small beakers, etc. Furnished with drop-in cover (part number 4403801).</p> |  <p>The diagram shows a rectangular stainless steel basket with a grid-like base and vertical side rails. A separate, matching grid-like cover is shown above the basket, designed to fit inside it.</p>                  |
| 4591601                | <p><b>48-Pin Insert – Stainless Steel.</b> For use with beakers, Erlenmeyer flasks, and wide-mouth glassware. For placement in bottom or upper rack.</p>                      |  <p>The diagram illustrates a rectangular stainless steel insert with a grid base. It features 48 vertical pins of varying heights, arranged in a regular pattern to hold glassware like beakers and flasks upright.</p> |
| 4587000<br><br>4595900 | <p><b>Top Rack – Stainless Steel.</b> Included with SteamScrubbers.</p> <p><b>Top Rack – Stainless Steel.</b> Includes 2 slides for installation in FlaskScrubber.</p>        |  <p>The diagram shows a rectangular stainless steel top rack with a grid base and vertical side rails. Two long, cylindrical slides are shown resting on the grid, used for installation into a flask scrubber.</p>     |
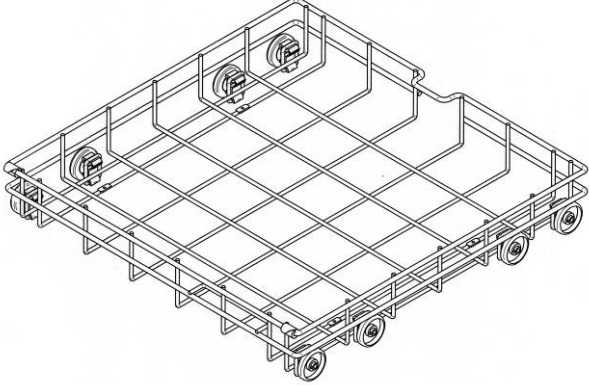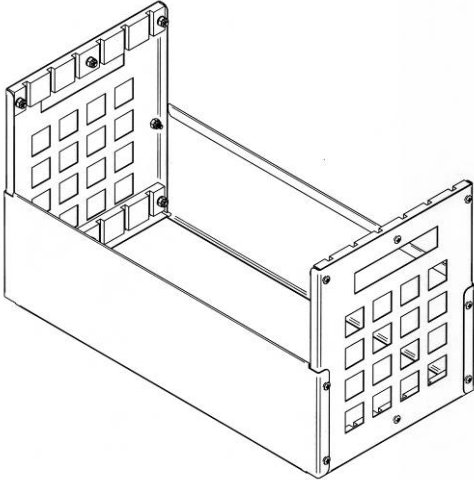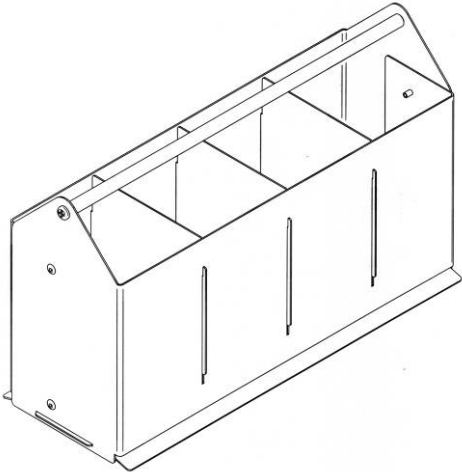
| PART #  | DESCRIPTION  |  |
|---------|--|--|
| 4588500 | <p><b>Bottom Rack – Stainless Steel.</b> Included with SteamScrubber.</p>  |    |
| 4591500 | <p><b>DNA Sequencing Plate Insert.</b> For electrophoresis plates 10.0 x 16.6 inches (25.4 x 42.1 cm). Holds 6 plates.</p> |   |
| 4542500 | <p><b>Utensil Holder.</b> For spatulas, spoons, scoops &amp; miscellaneous utensils.</p>                                   |  |
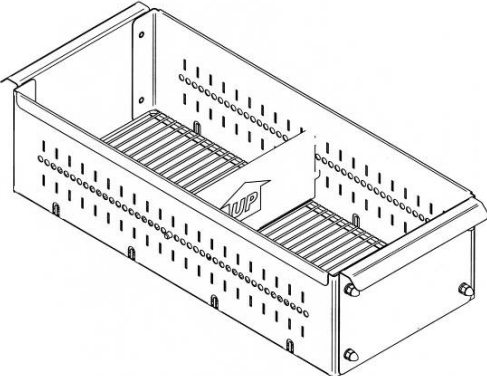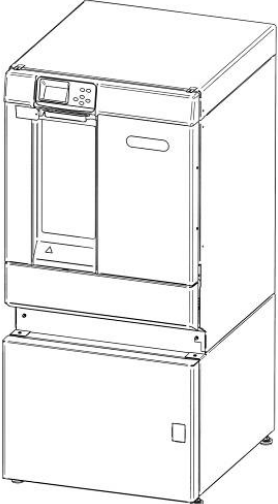
| PART #                                     | DESCRIPTION  |  |
|--|--|--|
| 4542100                                    | <b>Bulk Tube Insert.</b> For vertically positioning test tubes.  |    |
| 4595400                                    | <b>Stand.</b> For use with Freestanding models. Raises washer approximately 17 inches (43 cm). Front opens to provide storage underneath washer. Shown with glassware washer installed.        |  |
| 4591900<br>(115V)<br><br>4591901<br>(230V) | <b>Cool Drain Water Drain Kit.</b> Provides cool water to the drain plumbing. Connects to a cold water source.   |  |
| 4587500                                    | <b>Liquid Detergent Dispense Kit.</b> Permits liquid detergent to be dispensed into the glassware washer directly from a 1 gallon container.   |  |
| 4596100<br>(115V)<br><br>4596101<br>(230V) | <b>Mobile Conversion Kit.</b> Permits a Freestanding Glassware Washer to be converted to a Mobile model. Includes casters, electrical cord and plug, and plumbing to connect to a sink faucet. |  |

| <b>PART #</b> | <b>DESCRIPTION</b>   |  |
|---------------|--|--|
| 4592200       | <b>Connection Kit.</b> DI Water inlet valve to plastic tubing. |  |
| 1233700       | <b>Power Boost Transformer.</b> For 208V operation.            |  |


## **CONSUMABLES**

| <b>PART #</b> | <b>DESCRIPTION</b>                                    |
|---------------|---|
| 4422000       | 10-lb. (4.5 kg) LabSolutions Powder Detergent         |
| 4422100       | 27.5-lb. (12.5 kg) LabSolutions Powder Detergent      |
| 4522000       | 1 gallon (3.8 L) LabSolutions Liquid Detergent        |
| 4522200       | 1 liter (34 oz.) LabSolutions Neutralizing Acid Rinse |


# APPENDIX B

## GLASSWARE WASHER DIMENSIONS

### Freestanding Washer

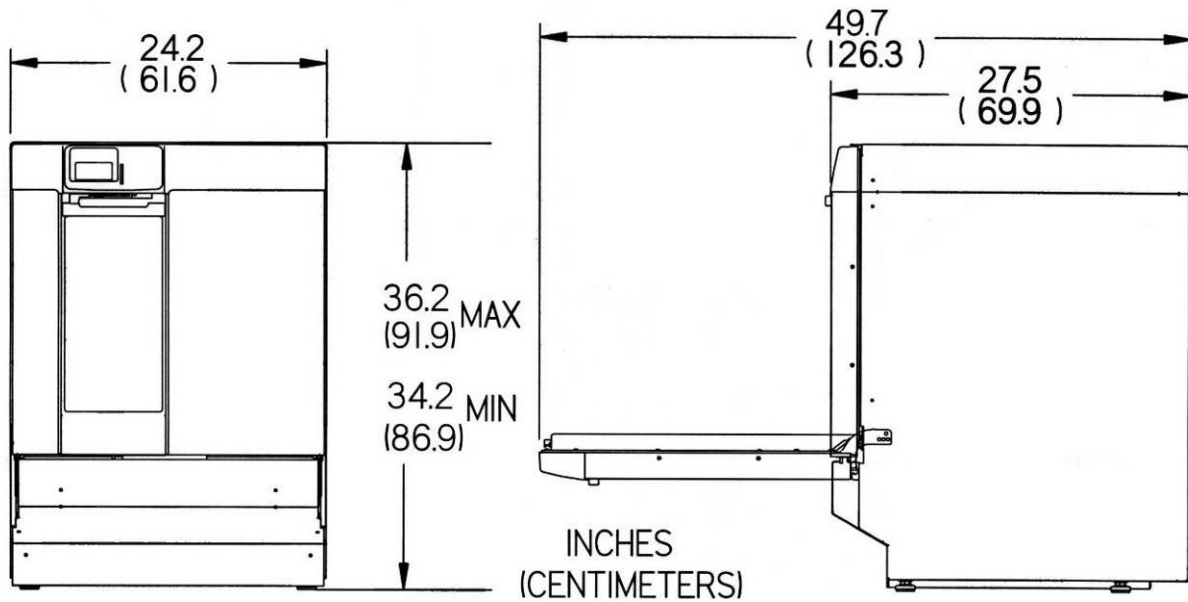


### Undercounter Washer

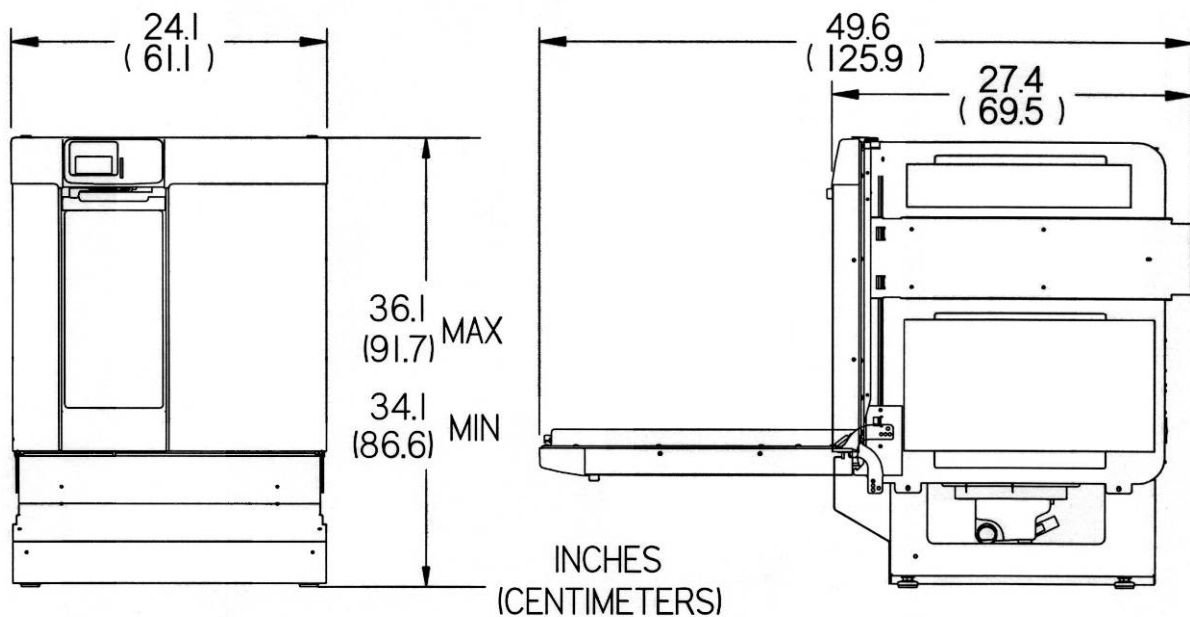




### Interior - All SteamScrubbers

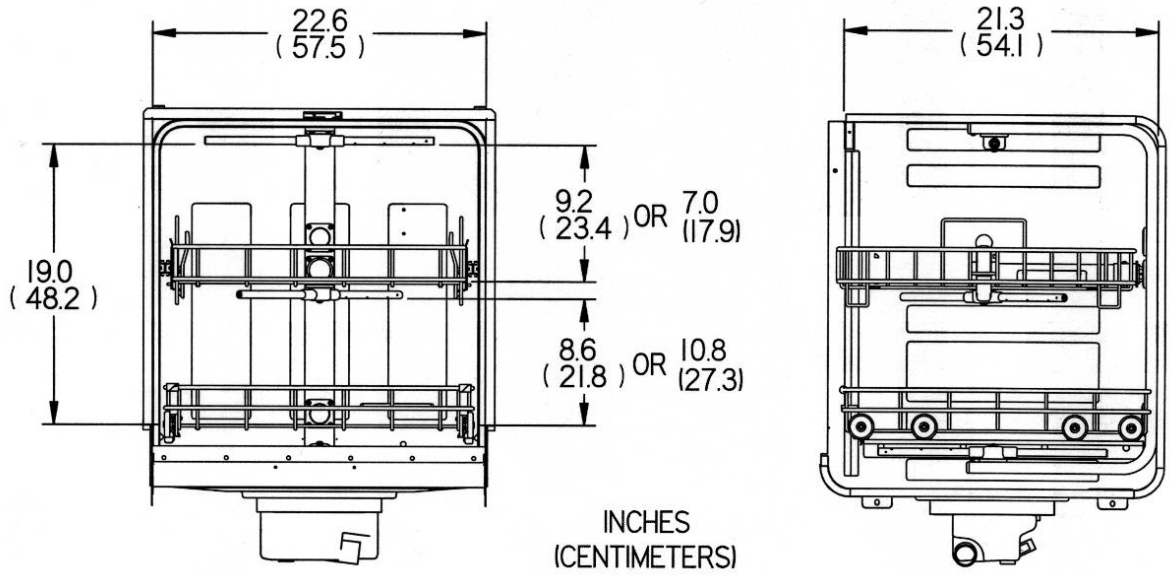

### Interior - All FlaskScrubbers

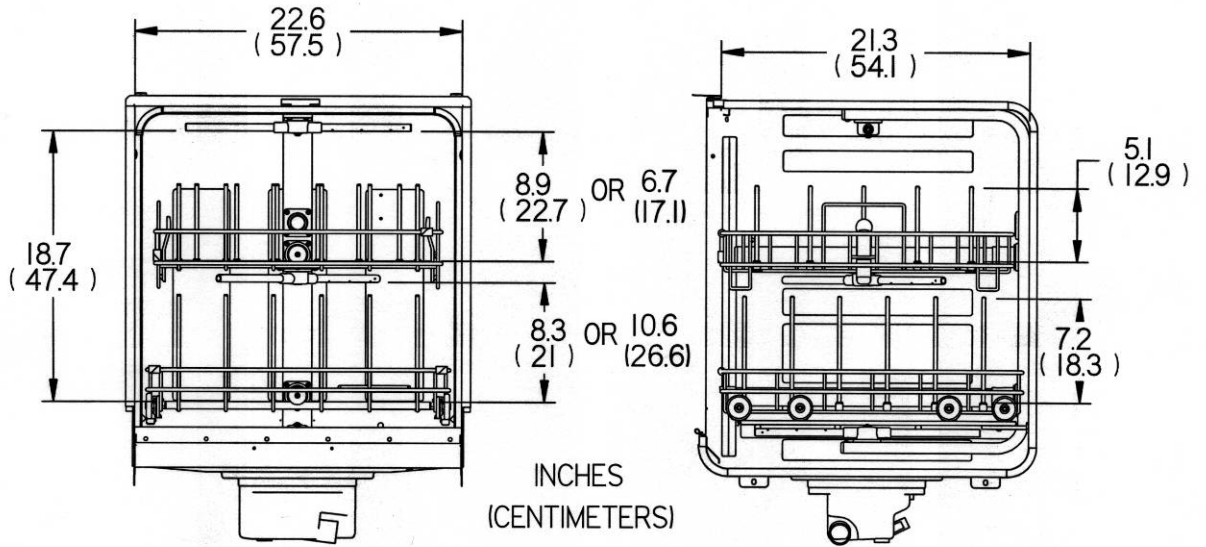

---

# APPENDIX C

## GLASSWARE WASHER

### SPECIFICATIONS

This Appendix contains various specifications for the Glassware Washer, including program times and water consumption for both the 115V and 230V models.

#### **Glassware Washer Specifications**

- Minimum fill water temperature: 120°F (49°C)
- Water consumption per fill: 3.4 gallons (13 liters)
- Minimum feed water pressure: 20 psi (138 kPa)
- Maximum feed water pressure: 120 psi (827 kPa)
- Minimum purified water feed pressure: 0 psi (0 kPa)
- Heat rejected from washer: less than 104 BTU/minute (115V models); less than 170 BTU/minute (230V models)
- Water heating rate: 1.6°F/minute (0.9°C/minute) 115V models operating on 115V; 2.2°F/minute (1.2°C/minute) 230V models operating on 208V

#### **Electrical Specifications**

- Operating voltage for 115V washers: 104/127V
- Operating voltage for 230V washers: 207/253V
- Maximum amperage for 115 volt washers: 16.0 amperes
- Maximum amperage for 230 volt washers: 12.0 amperes
- Frequency 115 volt washers: 60 Hz
- Frequency 230 volt washers: 50/60 Hz
- Phase: Single
- Sump element power: 1,200 watts (115V models); 2,000 watts (230V models)

### Environmental Conditions

- Indoor use only.
- Maximum altitude: 6562 feet (2000 meters).
- Ambient temperature range: 41° to 104°F (5° to 40°C).
- Maximum relative humidity: 80% for temperatures up to 88°F (31°C), decreasing linearly to 50% relative humidity at 104°F (40°C).
- Main supply voltage fluctuations not to exceed  $\pm 10\%$  of the nominal voltage.
- Transient overvoltages according to Installation Categories II (Overvoltage Categories per IEC 1010). Temporary voltage spikes on the AC input line that may be as high as 1500V for 115V models and 2500V for 230V models are allowed.
- Used in an environment of Pollution degrees 2 (i.e., where normally only non-conductive atmospheres are present). Occasionally, however, a temporary conductivity caused by condensation must be expected, in accordance with IEC 664.



Program Times & Water Consumption (115V Models)

Factory Program Descriptions and Possible User Modifications- 115V Models

|  | Units                 | <u>Rinse Only</u> |                             | <u>Plastic</u> |                              | <u>Glass</u> |                              | <u>Glass Plus</u> |                              | <u>Scientific</u> |                              |
|--|-----------------------|-------------------|-----------------------------|----------------|------------------------------|--------------|------------------------------|-------------------|------------------------------|-------------------|------------------------------|
|  |                       | Factory           | User set                    | Factory        | User set                     | Factory      | User set                     | Factory           | User set                     | Factory           | User set                     |
| Delayed Start  | (hours)               | 0                 | 1-8                         | 0              | 1-8                          | 0            | 1-8                          | 0                 | 1-8                          | 0                 | 1-8                          |
| Wash 1 Time  | (minutes)             |                   |                             | 0              | None                         | 5            | None                         | 5                 | None                         | 5                 | None                         |
| Steam  |                       |                   |                             | No             | None                         | No           | Yes / No                     | Yes               | Yes / No                     | No                | Yes / No                     |
| Wash 2 Time  | (minutes)             |                   |                             | 10             | 10,20,30,40                  | 10           | 10,20,30,40                  | 10                | 10,20,30,40                  | 20                | 10,20,30,40                  |
| Wash 2 Temp  | (°F)<br>(°C)          |                   |                             | 122<br>50      | 100 to 140<br>38-60          | 140<br>60    | 100 to 180<br>38-82          | 140<br>60         | 100 to 180<br>38-82          | 158<br>70         | 100 to 180<br>38-82          |
| Number of Rinses   |                       | 1                 | 0-6                         | 2              | 0-6                          | 3            | 0-6                          | 3                 | 0-6                          | 4                 | 0-6                          |
| No of Pure Rinses  |                       | 0                 | 0-6                         | 0              | 0-6                          | 0            | 0-6                          | 0                 | 0-6                          | 0                 | 0-6                          |
| Final Rinse  | (°F)<br>(°C)          | 100<br>38         | 100 to 140<br>38-60         | 122<br>50      | 100 to 140<br>38-60          | 140<br>60    | 100 to 180<br>38-82          | 140<br>60         | 100 to 180<br>38-82          | 158<br>70         | 100 to 180<br>38-82          |
| Dry Time   | (minutes)             | 0                 | 0 or 15-99                  | 15             | 0 or 15-99                   | 30           | 0 or 15-99                   | 30                | 0 or 15-99                   | 40                | 0 or 15-99                   |
| Dry Temp   | (°F)<br>(°C)          | 100<br>38         | 100-140<br>38-60            | 122<br>50      | 100 to 140<br>38-60          | 140<br>60    | 100 to 158<br>38-70          | 140<br>60         | 100 to 158<br>38-70          | 140<br>60         | 100 to 158<br>38-70          |
| Total Water Consumption  | (gallons)<br>(liters) | 3.4<br>12.9       | 3.4 to 20.4<br>12.9 to 77.2 | 13.6<br>51.5   | 6.8 to 27.2<br>25.7 to 103.0 | 17.0<br>64.4 | 6.8 to 27.2<br>25.7 to 103.0 | 17.0<br>64.4      | 6.8 to 27.2<br>25.7 to 103.0 | 20.4<br>77.2      | 6.8 to 27.2<br>25.7 to 103.0 |
| Total Wash Time<br>(Excluding delays for heating water and dry time) | (min)                 | 9                 |                             | 48             |                              | 81           |                              | 91                |                              | 110               |                              |

Note  
Times shown are for tap water rinses. Pure rinses add 1.5 minutes per rinse.  
Additional time is required to allow inlet water to heat up to the set temperature.  
Steam adds 10 minutes to the total program time.

**Factory Program Descriptions and Possible User Modifications- 115V Models**

|  | Units                 | <u>Scientific Plus</u> |                              |         | <u>Dry Only</u> |           |                              | <u>User 1</u> |                              |              | <u>User 2</u>                |         |          |
|--|-----------------------|------------------------|------------------------------|---------|-----------------|-----------|------------------------------|---------------|------------------------------|--------------|------------------------------|---------|----------|
|  |                       | Factory                | User set                     | Factory | User set        | Factory   | User set                     | Factory       | User set                     | Factory      | User set                     | Factory | User set |
| Delayed Start  | (hours)               | 0                      | 1-8                          |         |                 | 0         | 1-8                          | 0             | 1-8                          | 0            | 1-8                          |         |          |
| Wash 1 Time  | (minutes)             | 5                      | None                         |         |                 | 5         | None                         | 5             | None                         | 5            | None                         |         |          |
| Steam  |                       | Yes                    | Yes / No                     |         |                 | No        | Yes / No                     | No            | Yes / No                     | No           | Yes / No                     |         |          |
| Wash 2 Time  | (minutes)             | 20                     | 10,20,30,40                  |         |                 | 10        | 10,20,30,40                  | 10            | 10,20,30,40                  | 10           | 10,20,30,40                  |         |          |
| Wash 2 Temp  | (°F)<br>(°C)          | 158<br>70              | 100 to 180<br>38-82          |         |                 | 140<br>60 | 100 to 180<br>38-82          | 180<br>82     | 100 to 180<br>38-82          | 180<br>82    | 100 to 180<br>38-82          |         |          |
| Number of Rinses   |                       | 4                      | 0-6                          |         |                 | 2         | 0-6                          | 2             | 0-6                          | 2            | 0-6                          |         |          |
| No of Pure Rinses  |                       | 0                      | 0-6                          |         |                 | 0         | 0-6                          | 0             | 0-6                          | 0            | 0-6                          |         |          |
| Final Rinse  | (°F)<br>(°C)          | 158<br>70              | 100 to 180<br>38-82          |         |                 | 140<br>60 | 100 to 180<br>38-82          | 180<br>82     | 100 to 180<br>38-82          | 180<br>82    | 100 to 180<br>38-82          |         |          |
| Dry Time   | (minutes)             | 40                     | 0 or 15-99                   |         |                 | 60        | 0 or 15-99                   | 15            | 0 or 15-99                   | 15           | 0 or 15-99                   |         |          |
| Dry Temp   | (°F)<br>(°C)          | 140<br>60              | 100 to 158<br>38-70          |         |                 | 140<br>60 | 100 to 158<br>38-70          | 140<br>60     | 100 to 158<br>38-70          | 140<br>60    | 100 to 158<br>38-70          |         |          |
| Total Water Consumption  | (gallons)<br>(liters) | 20.4<br>77.2           | 6.8 to 27.2<br>25.7 to 103.0 |         |                 | 0<br>0    | 6.8 to 27.2<br>25.7 to 103.0 | 13.6<br>51.5  | 6.8 to 27.2<br>25.7 to 103.0 | 13.6<br>51.5 | 6.8 to 27.2<br>25.7 to 103.0 |         |          |
| Total Wash Time<br>(Excluding delays for heating water and dry time) | (min)                 | 120                    |                              |         |                 | 60        |                              | 58            |                              | 58           |                              |         |          |

Program Times & Water Consumption (230V Models)

Factory Program Descriptions and Possible User Modifications- 230V Models

|  | Units                 | Rinse Only  |                             | Plastic      |                              | Glass        |                              | Glass Plus   |                              | Scientific   |                              | Scientific Plus |                              |
|--|-----------------------|-------------|-----------------------------|--------------|------------------------------|--------------|------------------------------|--------------|------------------------------|--------------|------------------------------|-----------------|------------------------------|
|  |                       | Factory     | User set                    | Factory      | User set                     | Factory      | User set                     | Factory      | User set                     | Factory      | User set                     | Factory         | User set                     |
| Delayed Start  | (hours)               |             |                             | 0            | 0-8 Hr                       | 0            | 0-8 Hr                       | 0            | 0-8 Hr                       | 0            | 0-8 Hr                       | 0               | 0-8 Hr                       |
| Wash 1 Time  | (minutes)             |             |                             | 5            | None                         | 5            | None                         | 5            | None                         | 5            | None                         | 5               | None                         |
| Steam  |                       |             |                             | N            | None                         | N            | Y N                          | Y            | Y N                          | N            | Y N                          | Y               | Y N                          |
| Wash 2 Time  | (minutes)             |             |                             | 10           | 10,20,30,40                  | 10           | 10,20,30,40                  | 10           | 10,20,30,40                  | 20           | 10,20,30,40                  | 20              | 10,20,30,40                  |
| Wash 2 Temperature   | (°F)                  |             |                             | 122          | 100 to 140                   | 140          | 100 to 199                   | 140          | 100 to 199                   | 158          | 100 to 199                   | 158             | 100 to 199                   |
|  | (°C)                  |             |                             | 50           | 38-60                        | 60           | 38-93                        | 60           | 38-93                        | 70           | 38-93                        | 70              | 38-93                        |
| Number of Rinses   |                       | 1           | 0-6                         | 2            | 0-6                          | 3            | 0-6                          | 3            | 0-6                          | 4            | 0-6                          | 4               | 0-6                          |
| No of Pure Rinses  |                       | 0           | 0-6                         | 0            | 0-6                          | 0            | 0-6                          | 0            | 0-6                          | 0            | 0-6                          | 0               | 0-6                          |
| Final Rinse  | (°F)                  | 100         | 100 to 140                  | 122          | 100 to 140                   | 140          | 100 to 199                   | 140          | 100 to 199                   | 158          | 100 to 199                   | 158             | 100 to 199                   |
|  | (°C)                  | 38          | 38-60                       | 50           | 38-60                        | 60           | 38-93                        | 60           | 38-93                        | 70           | 38-93                        | 70              | 38-93                        |
| Dry Time   | (minutes)             | 0           | 0 or 15-99                  | 15           | 0 or 15-99                   | 30           | 0 or 15-99                   | 30           | 0 or 15-99                   | 40           | 0 or 15-99                   | 40              | 0 or 15-99                   |
| Dry Temperature  | (°F)                  | 100         | 100 to 140                  | 122          | 100 to 140                   | 140          | 100 to 158                   | 140          | 100 to 158                   | 140          | 100 to 158                   | 140             | 100 to 158                   |
|  | (°C)                  | 38          | 38-60                       | 50           | 38-60                        | 60           | 38-70                        | 60           | 38-70                        | 60           | 38-70                        | 60              | 38-70                        |
| Total Water Consumption  | (gallons)<br>(liters) | 3.4<br>12.9 | 3.4 to 20.4<br>12.9 to 77.2 | 13.6<br>51.5 | 6.8 to 27.2<br>25.7 to 103.0 | 17.0<br>64.4 | 6.8 to 27.2<br>25.7 to 103.0 | 17.0<br>64.4 | 6.8 to 27.2<br>25.7 to 103.0 | 20.4<br>77.2 | 6.8 to 27.2<br>25.7 to 103.0 | 20.4<br>77.2    | 6.8 to 27.2<br>25.7 to 103.0 |
| Total Wash Time<br>(Excluding delays for heating water and dry time) | (min)                 | 9           |                             | 48           |                              | 81           |                              | 91           |                              | 110          |                              | 120             |                              |

Note

Times shown are for tap water rinses. Pure rinses add 1.5 minutes per rinse. Additional time is required to allow inlet water to heat up to the set temperature. Steam adds 10 minutes to the total program time.

**Factory Program Descriptions and Possible User Modifications- 230V Models**

|  | Units     | <u>Intense</u> |               | <u>Intense Plus</u> |               | <u>Extreme</u> |               | <u>Dry Only</u> |             | <u>User 1</u> |               | <u>User 2</u> |               |
|--|-----------|----------------|---------------|---------------------|---------------|----------------|---------------|-----------------|-------------|---------------|---------------|---------------|---------------|
|  |           | Factory        | User set      | Factory             | User set      | Factory        | User set      | Factory         | User set    | Factory       | User set      | Factory       | User set      |
| Delayed Start  | (hours)   | 0              | 0-8 Hr        | 0                   | 0-8 Hr        | 0              | 0-8 Hr        | 0               | 0-8 Hr      | 0             | 0-8 Hr        | 0             | 0-8 Hr        |
| Wash 1 Time  | (minutes) | 10             | None          | 10                  | None          | 10             | None          | 5               | None        | 5             | None          | 5             | None          |
| Steam  |           | N              | Y N           | Y                   | Y N           | Y              | Y N           | N               | Y N         | N             | Y N           | N             | Y N           |
| Wash 2 Time  | (minutes) | 30             | 10,20,30,40   | 30                  | 10,20,30,40   | 40             | 10,20,30,40   | 10              | 10,20,30,40 | 10            | 10,20,30,40   | 10            | 10,20,30,40   |
| Wash 2 Temperature   | (°F)      | 180            | 100 to 199    | 180                 | 100 to 199    | 199            | 100 to 199    | 140             | 100 to 199  | 140           | 100 to 199    | 180           | 100 to 199    |
|  | (°C)      | 82             | 38-93         | 82                  | 38-93         | 93             | 38-93         | 60              | 38-93       | 60            | 38-93         | 82            | 38-93         |
| Number of Rinses   |           | 5              | 0-6           | 5                   | 0-6           | 6              | 0-6           | 2               | 0-6         | 2             | 0-6           | 2             | 0-6           |
| No of Pure Rinses  |           | 0              | 0-6           | 0                   | 0-6           | 0              | 0-6           | 0               | 0-6         | 0             | 0-6           | 0             | 0-6           |
| Final Rinse  | (°F)      | 180            | 100 to 199    | 180                 | 100 to 199    | 199            | 100 to 199    | 140             | 100 to 199  | 140           | 100 to 199    | 180           | 100 to 199    |
|  | (°C)      | 82             | 38-93         | 82                  | 38-93         | 93             | 38-93         | 60              | 38-93       | 60            | 38-93         | 82            | 38-93         |
| Dry Time   | (minutes) | 50             | 0 or 15-99    | 50                  | 0 or 15-99    | 60             | 0 or 15-99    | 60              | 0 or 15-99  | 15            | 0 or 15-99    | 15            | 0 or 15-99    |
| Dry Temperature  | (°F)      | 140            | 100 to 158    | 140                 | 100 to 158    | 140            | 100 to 158    | 140             | 100 to 158  | 140           | 100 to 158    | 140           | 100 to 158    |
|  | (°C)      | 60             | 38-70         | 60                  | 38-70         | 60             | 38-70         | 60              | 38-70       | 60            | 38-70         | 60            | 38-70         |
| Total Water Consumption  | (gallons) | 23.8           | 6.8 to 27.2   | 23.8                | 6.8 to 27.2   | 27.2           | 6.8 to 27.2   | 0               | 0           | 13.6          | 6.8 to 27.2   | 13.6          | 6.8 to 27.2   |
|  | (liters)  | 90.1           | 25.7 to 103.0 | 90.1                | 25.7 to 103.0 | 103.0          | 25.7 to 103.0 | 0               | 0           | 51.5          | 25.7 to 103.0 | 51.5          | 25.7 to 103.0 |
| Total Wash Time<br>(Excluding delays for heating water and dry time) | (min)     | 143            |               | 153                 |               | 182            |               | 60              |             | 58            |               | 58            |               |



**DECLARATION OF CONFORMITY**

Application Council Directive(s): 73/23/EEC, 89/336/EEC, 2002/95/EC (ROHS),  
2002/96/EC (WEEE), 2004/108/EC

Standard(s) to which conformity is declared: EN61010-1, EN61326-1, EN55022,  
EN61000-3-2/3

Manufacturer’s Name: Labconco Corporation

Manufacturer’s Address: 8811 Prospect Avenue  
Kansas City, MO 64132 USA

Importer’s Name: See Shipping/Customs Documents

Importer’s Address: See Shipping/Customs Documents for your equipment

Type of Equipment: Laboratory Equipment – Glassware Washers

Model No.: Washers with Micro Processor Control  
4400321 – 208-240V, 50/60 Hz – SteamScrubber Undercounter  
4400331 – 208-240V, 50/60 Hz – SteamScrubber Undercounter with glass window  
4400421 – 208-240V, 50/60 Hz – SteamScrubber Freestanding  
4400431 – 208-240V, 50/60 Hz – SteamScrubber Freestanding with glass window  
4400461 – 208-240V, 50/60 Hz – SteamScrubber Freestanding with electronics cleaning options  
4420321 – 208-240V, 50/60 Hz – FlaskScrubber Undercounter  
4420331 – 208-240V, 50/60 Hz – FlaskScrubber Undercounter with glass window  
4420421 – 208-240V, 50/60 Hz – FlaskScrubber Freestanding  
4420431 – 230V, 50/60 Hz – FlaskScrubber Freestanding with glass window  
4540031 – 208-240V, 50/60 Hz – FlaskScrubber Vantage (has built-in water conductivity probe)  
with glass window

Serial No.: Various – See Individual Declaration

Year of Manufacture: 2007 and Subsequent

I, the undersigned, hereby declare that the equipment specified above conforms to the  
above Directive(s) and Standard(s).

See individual Declaration of Conformity which  
will be signed by the importer for your country.

Place: \_\_\_\_\_  
(Signature)

Date: \_\_\_\_\_  
(Full Name)

\_\_\_\_\_  
(Position)

Labconco P/N 36960-61, REV. A, ECO E402